

Purification of Peptides by Twin-column Countercurrent Chromatography

Thomas Müller-Späth^{1,2}

¹Institute for Chemical and Bioengineering, ETH Zurich, 8093 Zurich, Switzerland

²ChromaCon AG, 8005, Zurich, Switzerland

TIDES conference, Amsterdam, 8th Nov 2018

Presentation Highlights

MCSGP is a counter-current chromatography process that simultaneously achieves high yield and purity in difficult peptide purifications.

- The increased yield of MCSGP ...
 - ✓ allows downscaling of the upstream chemical synthesis steps
 - improves productivity of the downstream process leading to smaller columns required
 - ✓ reduces solvent consumption
 - ✓ eliminates the need for re-chromatography
 - ✓ avoids generation of side-fractions to be stored and analyzed (reduction of analytical burden)

Presentation Outline

- Multicolumn chromatography (MCSGP) process introduction
- MControl Dynamic process control for MCSGP
- MCSGP for peptide purification case studies
- Economic evaluation of MCSGP



MCSGP -<u>Multicolumn Countercurrent Solvent Gradient</u> <u>Chromatography</u>

© ChromaCon AG 2018

Purification Challenge

Yield-purity trade-off for ternary separations



Background – MCSGP process principle

- Single column batch chromatography suffer from a yield-purity trade-off due to overlapping impurities:
- High purity can only be obtained at the cost of yield and vice versa
- Trade-off becomes worse with increasing load and increasing flow rate
- → In batch chromatography: Conflicting aims: purity vs. yield, load, productivity
- MCSGP can obtain high purity and yield simultaneously



Background – MCSGP process principle

- MCSGP (Multicolumn counter-current solvent gradient purification) is a chromatography process that uses two columns of the same type
- MCSGP uses internal recycling and inline dilution fractions to automatically recover the product from impure side-fractions, eluting only product of high purity at a high yield



MCSGP explained







MCSGP explained



MCSGP explained











MCSGP explained





Start over



MControl - Dynamic process control for MCSGP

© ChromaCon AG 2018

MControl – Robust MCSGP operation

In MCSGP, operation with MControl significantly reduces effects of the following parameters on product quality:

- temperature
- solvent quality
- conductivity, pH
- column variability (bed height, aging, packing quality)

MControl compensates for peak shifts by adjusting the fractionation start:

- →Same product fraction position
- → Same product quality
- → Increased robustness of continuous process operation



20

| 21

MCSGP adjusts fraction collection

MControl: Dynamic MCSGP Process Control

- Example: MCSGP run on Contichrom with two different columns
- Chromatograms show 6 cycles superimposed, small protein model system, cation-exchange, linear gradient elution
 - MControl runs the linear gradient at the same slope, prolonging the elution (t1, t2) until the UV threshold is reached and product collection starts (P).



Institute for Chemical and Bioengineering Morbidelli Group

MCSGP operation

- Cyclic steady state of MCSGP: Constant product concentration and purity
- MControl supports robust operation



Morbidelli Group

MCSGP for peptide purification – case studies

Peptide Case study*



Batch reference run



Gradient elution on Kromasil C-18. 10 um, 0.46 x 25 cm, solvents: water, ACN

Institute for Chemical and Bioengineering Morbidelli Group

Peptide Case study results

MCSGP resolved the yield/purity trade-off problem of batch:

- ✓ 70% yield increase at target purity
- ✓ 10x productivity improvement
- ✓ 70% decrease in solvent consumption (S.C.) at target purity



ChromaCon

Bristol-Myers Squibb

K A I

Case: Therapeutic peptide purification by MCSGP (Liraglutide)

LUND UNIVERSITY LUND UNIVERSITY LIBRARIES	Register publications S	t
Home Publications Departments Simulation, optimization and implementat of Liraglutide	new search	novo nordisk [®]
Gomis Fons, Joaquín ^{Lu} (2017) KET920 20171 Chemical Engineering	Mark	

- Collaboration between University of Lund (Sweden) and Novo Nordisk
- Detailed results are confidential
- Outcome: At high purity (98%, 99%), MCSGP was more favorable than batch chromatography
- Productivity was improved through MCSGP

Economic evaluation of MCSGP



Assumptions for economic evaluation

- Batch chromatography has varying yields of 40, 50, 60, 70% representing varying impurity content / purifications of peptides with different sizes (15, 20, 25-mer)
- Rationale: shorter peptides → less complex synthesis → fewer impurities → higher yield in batch chromatography
- Yield of MCSGP is 95%, independent of difficulty of separation / peptide size
- The load for all batch runs and MCSGP was assumed to be 10 g crude/L of stationary phase



Institute for Chemical and Bioengineering Morbidelli Group

Further Assumptions

Parameters for simulation of production and purification of 10 kg peptide:

		Batch	MCSGP
Column bed height	[cm]	25.0	10.0
Replacement of stationary phase	[%/year]	30	30
Stat. phase costs	[US\$/kg]	7000	7000
Synthesis batch size	[kg]	1	1
Synthesis costs / g	[US\$/g]	200	200
Synthesis costs / batch	[US\$]	200,000	200,000
Solvent costs	[US\$/L]	6	6
Chrom. system costs	[US\$]	500,000	1,500,000
Depreciation period	[a]	10	10
number of samples to be analyzed per cycle	[-]	10	1
QA/QC costs per sample	[US\$]	200	200
Plant operating costs	[US\$/day]	8,000	5,000
Max. time permitted for chromatography	[hrs]	16	16

 Further assumption for batch chromatography: Use of re-chromatography: 25% of yield loss can be recovered

Further assumptions

 Parameters for simulation of production and purification of 10 kg peptide:

Parameter	Unit	Batch 1	Batch 2	Batch 3	Batch 4	MCSGP
Yield	[%]	40	50	60	70	95
Flow rate	[cm/h]	181	181	181	181	271
Cycle time	[min]	232.7	232.7	232.7	232.7	80

- Assumed linear flow rate in MCSGP 50% higher than batch flow rate:
 - MCSGP can achieve high product yield in spite of a high flow rate which causes larger overlaps of product and impurities, due to its internal recycling capabilities.
 - the shorter bed height of MCSGP allows larger linear flow rates due to reduced backpressure.

Results: Total costs including synthesis and re-chromatography

Cost benefit of MCSGP: US\$ 0.6 million to US\$ 2.1 million / 10 kg peptide



Results: Cost difference batch - MCSGP

- Due to low yield of batch chromatography, additional synthesis batches need to be produced, driving up costs compared to MCSGP.
- With increasing chromatography yield the number of required extra batches decreases, improving overall costs



Results: Payback period of MCSGP

- Payback period of MCSGP is calculated relative to the batch runs with 40%, 50%, 60% and 70% yield respectively (10 kg of peptide per year)
- Payback = (CAPEX difference batch-MCGSP)/(savings through MCSGP per year)
- → In all cases the payback period of MCSGP is equal or less than 19 months
- → Payback period can be even shorter if more than 10 kg peptide is produced per year



Morbidelli Group

Results: Total costs including synthesis and re-chromatography

- Only chromatography cost shown (10 kg peptide):
 - Solvent cost dominating for batch chromatography
 - CAPEX dominating for MCSGP
- Savings through MCSGP (chromatography costs only) : US\$ 200k US\$ 500k



Results: Solvent consumption

- MCSGP cuts solvent consumption by up to 85%, corresponding to 56,000 L p.a.
- Additional cost savings through reduced solvent handling/storage/disposal /recycling ... not included in evaluation but in favor of MCSGP



Results: Column and pump sizes

Comparison of Batch and MCSGP:

		Batch 1	Batch 2	Batch 3	Batch 4	MCSGP
Yield	[%]	40	50	60	70	95
Column inner diameter	[cm]	60	60	60	60	30
Column volume	[L]	70.7	70.7	70.7	70.7	2x 7.1
Required pump size on skid	[L/min]	8.5	8.5	8.5	8.5	3.2

- Column diameter reduced from 60 cm i.d. to 2x 30 cm i.d.
- Total column packing volume reduced from 70.7 L to 14.2 L
- Pump flow rate on skid reduced from 8.5 L/min (510 L/h) to 3.2 L/min (190 L/h)



Results: Chromatography costs

- Sensitivity analysis: use of smaller columns.
- Use of smaller columns reduces stationary phase costs but drives up QA/QC costs and plant operating cost through the increased processing time → increased chromatography costs



Morbidelli Group

MCSGP scalability

Lab Scale systems

Cost-competitive, all-in-one process capabilities

Pilot / Production-scale (GMP)

• High throughput, reduced costs



Contichrom CUBE 30/100



Contichrom TWIN MCSGP HPLC GMP systems





Institute for Chemical and Bioengineering Morbidelli Group Flow rates: 100 mL/min

Flow rates: up to 20 L/min

Conclusions

- MCSGP simultaneously achieves high yield and purity in difficult peptide purifications
- The increased yield of MCSGP ...
 - ✓ allows downscaling of the upstream chemical synthesis steps
 - ✓ improves productivity of the downstream process leading to smaller columns required
 - ✓ reduces solvent consumption
 - ✓ eliminates the need for re-chromatography
 - avoids generation of side-fractions to be stored and analyzed (reduction of analytical burden)
- All abovementioned points lead to massive cost savings compared to single column batch chromatography
- Economic analysis: Savings for an annual production amount of 10 kg peptide (synthesis, chromatography and re-chromatography) from US\$ 0.6 million to US\$ 2.1 million expected, in comparison to the single column reference process

Thank you for attending any questions?

