

# YMC-Triart Bio C4

YMC-Triart Bio C4 is a novel, organic/inorganic hybrid silica derivatized with C4 (L26 USP designation) based on wide-pore (300Å) particles. This stationary phase is designed for HPLC and UHPLC separations of proteins and oligonucleotides, and is particularly useful for intact antibody analyses run at high temperatures.

Triart Bio C4 is well-suited for high-sensitivity analysis using mass spec-friendly mobile phases containing formic acid and/or TFA. As with all Triart products, chromatographers enjoy wide temperature and pH ranges (up to 90°C at pH 1-7, and 50°C at pH 7-10), allowing these parameters to be used as tools for optimizing separation conditions without having to worry about stationary phase durability.


## Features

- Designed for the separation of intact antibodies
- Available in 3 and 5µm for HPLC, and 1.9µm for UHPLC
- Stable over a wide pH range (1-10) and at high temperatures (up to 90°C)
- Proprietary end-capping makes for excellent peak shapes of proteins
- The chemical and physical stability of the Triart base material contributes to long column lifetime
- Rigorous manufacturing controls result in legendary YMC column-to-column and lot-to-lot reproducibility

### Specifications

Base	: Organic/inorganic hybrid silica
Stationary Phase	: C4 (USP L26)
Particle size	: 1.9, 3, 5µm
Pore size	: 300Å
Bonding	: Polymeric type
Endcapping	: Yes
Usable pH range	: 1-10
Max. temperature	: 90°C (pH 1-7) 50°C (pH 7-10)

## Superior performance for LC/MS



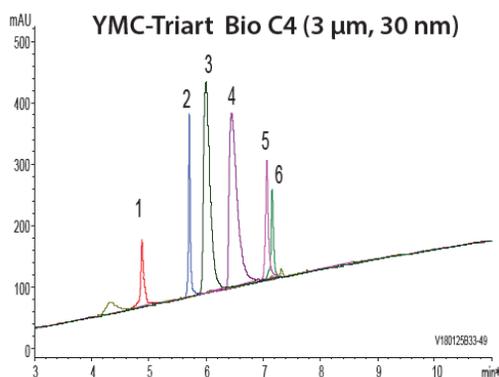



Sample:

1. Cytochrome-C (Horse heart)
2. Insulin (Bovine pancreas)
3. Transferrin (Human)
4. BSA
5. β-Lactoglobulin (Bovine)
6. α-Chymotrypsinogen A (Bovine pancreas)

Eluent	: A) water/formic acid (100/0.1) B) acetonitrile/formic acid (100/0.1) 10-95%B (0-15 min)
Temperature	: 40°C
Detection	: UV at 220nm

Triart Bio C4 exhibits improved peak shape and recovery with mobile phases containing formic acid/TFA, which are commonly used in LC/MS methods. This makes Triart Bio C4 ideal for high-sensitivity analysis of proteins.

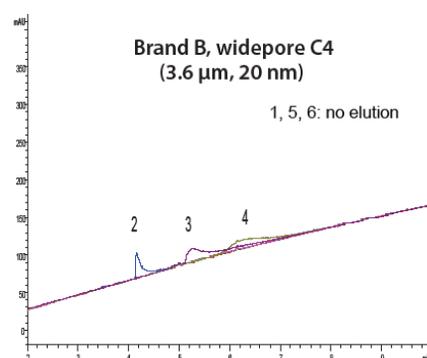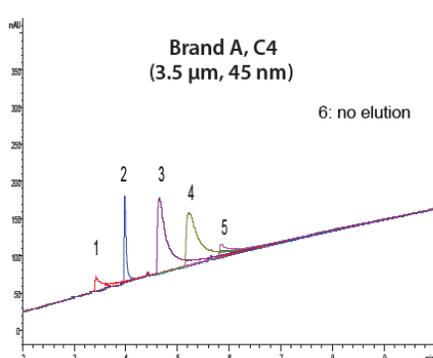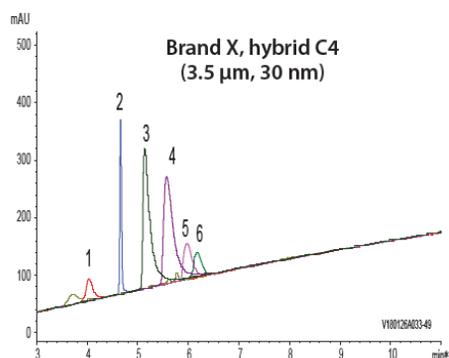

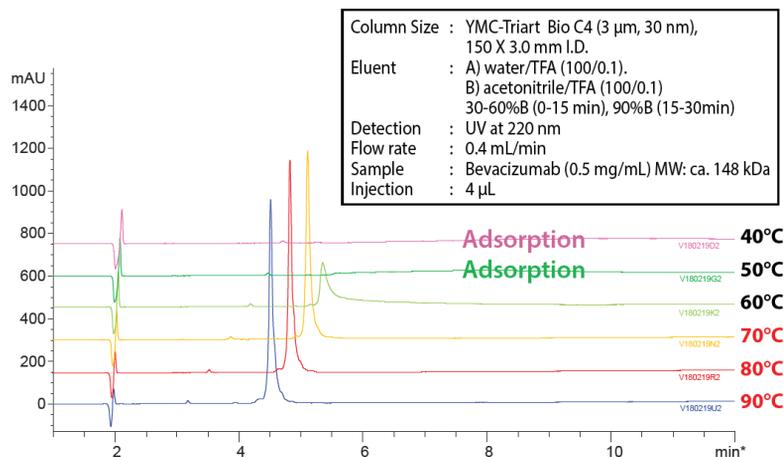

## Separation of Bevacizumab

When analyzing monoclonal antibodies by reversed phase, high temperature analysis is often required to minimize adsorption and obtain good peak shape. YMC-Triart Bio C4 with its high temperature and chemical durability is ideal for use at these conditions.

Improved peak shape was achieved at 60°C and higher, whereas no elution occurred at 50°C and below.

## Ordering Information

Particle Size (μm)	Column I.D. x length (mm)	Stainless Steel Column Product Number	PEEK-Lined Column Product Number
1.9	2.1 x 20	TB30SP9-02Q1PT	
	2.1 x 30	TB30SP9-03Q1PT	
	2.1 x 50	TB30SP9-05Q1PT	TB30SP9-05Q1PTP
	2.1 x 75	TB30SP9-L5Q1PT	
	2.1 x 100	TB30SP9-10Q1PT	TB30SP9-10Q1PTP
	2.1 x 150	TB30SP9-15Q1PT	TB30SP9-15Q1PTP
	3.0 x 50	TB30SP9-0503PT	
	3.0 x 75	TB30SP9-L503PT	
	3.0 x 100	TB30SP9-1003PT	
3.0	2.1 x 20	TB30S03-02Q1PTH	
	2.1 x 33	TB30S03-H3Q1PTH	
	2.1 x 50	TB30S03-05Q1PTH	TB30S03-05Q1PTP
	2.1 x 75	TB30S03-L5Q1PTH	
	2.1 x 100	TB30S03-10Q1PTH	TB30S03-10Q1PTP
	2.1 x 150	TB30S03-15Q1PTH	TB30S03-15Q1PTP
	3.0 x 50	TB30S03-0503PTH	
	3.0 x 75	TB30S03-L503PTH	
	3.0 x 100	TB30S03-1003PTH	
	3.0 x 150	TB30S03-1503PTH	
	4.6 x 33	TB30S03-H346PTH	
	4.6 x 50	TB30S03-0546PTH	TB30S03-0546PTP
	4.6 x 75	TB30S03-L546PTH	
	4.6 x 100	TB30S03-1046PTH	TB30S03-1046PTP
	4.6 x 150	TB30S03-1546PTH	TB30S03-1546PTP
	4.6 x 250	TB30S03-2546PTH	

Particle Size (μm)	Column I.D. x length (mm)	Stainless Steel Column Product Number	PEEK-Lined Column Product Number
5.0	2.1 x 20	TB30S05-02Q1PTH	
	2.1 x 33	TB30S05-H3Q1PTH	
	2.1 x 50	TB30S05-05Q1PTH	TB30S05-05Q1PTP
	2.1 x 75	TB30S05-L5Q1PTH	
	2.1 x 100	TB30S05-10Q1PTH	TB30S05-10Q1PTP
	2.1 x 150	TB30S05-15Q1PTH	TB30S05-15Q1PTP
	3.0 x 50	TB30S05-0503PTH	
	3.0 x 75	TB30S05-L503PTH	
	3.0 x 100	TB30S05-1003PTH	
	3.0 x 150	TB30S05-1503PTH	
	4.0 x 150	TB30S05-1504PTH	
	4.0 x 250	TB30S05-2504PTH	
	4.6 x 33	TB30S05-H346PTH	
	4.6 x 50	TB30S05-0546PTH	TB30S05-0546PTP
	4.6 x 75	TB30S05-L546PTH	
	4.6 x 100	TB30S05-1046PTH	TB30S05-1046PTP
4.6 x 150	TB30S05-1546PTH	TB30S05-1546PTP	
4.6 x 250	TB30S05-2546PTH		

Inquire about other column sizes.

Maximum pressure: 100MPa for 1.9μm;  
45 MPa for 3 and 5μm

Endfitting: Parker style (UHPLC compatible)

### Worldwide Availability

YMC America, Inc.  
www.ymcamerica.com

YMC Europe GmbH  
www.ymc.de

YMC Switzerland LLC  
www.ymc-schweiz.ch

YMC Shanghai Rep. Office  
www.ymcchina.com

YMC India Pvt. Ltd.  
www.ymcindia.com

YMC Korea Co., Ltd.  
www.ymckorea.com

YMC Taiwan Co., Ltd.  
www.ymctaiwan.com

YMC Singapore Tradelinks Pte. Ltd.  
www.ymc.co.jp/en/



## YMC CO., LTD.

YMC Karasuma-Gojo Bldg., 284 Daigo-cho  
Karasuma Nishiiru Gojo-dori, Shimogyo-ku  
Kyoto, 600-8106, Japan  
TEL:+81-75-342-4515 FAX:+81-75-342-4550  
www.ymc.co.jp sales@ymc.co.jp

Distributor

## YMC America, Inc.

8 Charlestown Street  
Devens, MA 01434  
Tel: +1 978 487 1100  
info@ymcamerica.com  
www.ymcamerica.com