


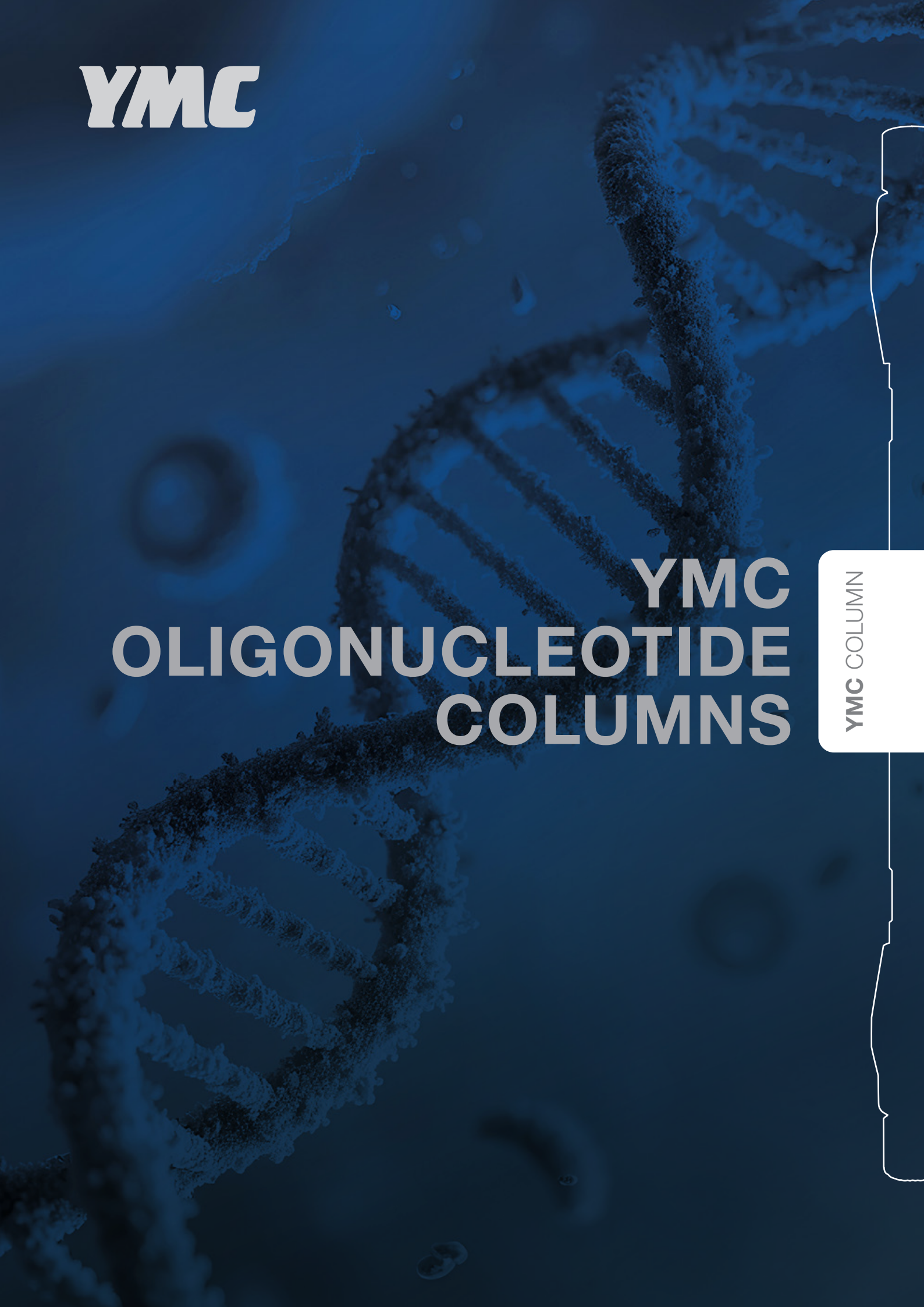

YMC

**YMC
OLIGONUCLEOTIDE
COLUMNS**

YMC COLUMN

YMC OLIGONUCLEOTIDE COLUMNS

YMC COLUMN



YMC



CONTENT

PAGE

Oligonucleotide Applications	07-12
IP-RP	13-18
Accura Bioinert Hardware	19-24
AEX	25-28
SEC	29-32
HILIC	33-36
Ordering Information	37-42

NUCLEIC ACIDS

DIFFERENT TYPES OF THERAPEUTIC NUCLEIC ACIDS

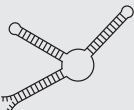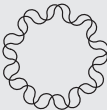	Antisense oligonucleotide/ miRNA	siRNA	Aptamer	gRNA	mRNA	Plasmid
Example structure						
Type(s) of nucleic acids	ssDNA, ssRNA	dsRNA	ss/dsRNA, ss/dsDNA	ss/dsRNA	ssRNA	dsDNA
Length (mer)	11–24	19–23	20–100	100	1,000–7,000	4,000–10,000
Mass range (kDa)	5	15	5–30	30	300–2,000	2,000–7,000
Modifications (phosphates, ribose, nucleobase)	Yes	Yes	Yes	Yes	Nucleobases only (5' cap, 3' Poly(A) tail)	No
Chemical conjugates / carrier systems	Yes / carrier possible	Yes / carrier possible	Yes / carrier possible	No / carrier possible	No / Lipid nanoparticles	No / carrier possible
Mode of action	mRNA degradation	mRNA degradation	Protein function modulation	Genome editing	Protein production	Gene expression

NUCLEIC ACIDS





PHASE SELECTION GUIDE

OLIGONUCLEOTIDES NUCLEIC ACIDS			
IP-RP	AEX	SEC	HILIC
YMC-Triart C18	BioPro IEX QF	YMC-Pack Diol-120	YMC-Triart Diol-HILIC
YMC-Triart Bio C18		YMC-Pack Diol-200	
Hydrosphere C18		YMC-Pack Diol-300	
YMC-Triart C8			
YMC-Triart Bio C4			




YMC OLIGONUCLEOTIDE APPLICATIONS

YMC COLUMN

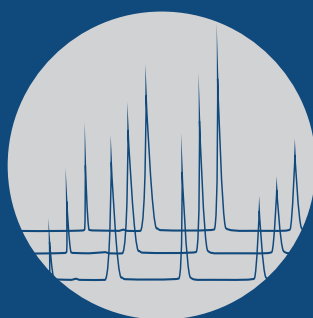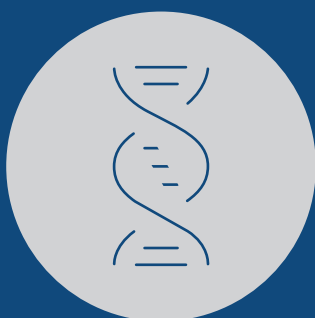

OLIGONUCLEOTIDE APPLICATIONS

Synthetic oligonucleotides

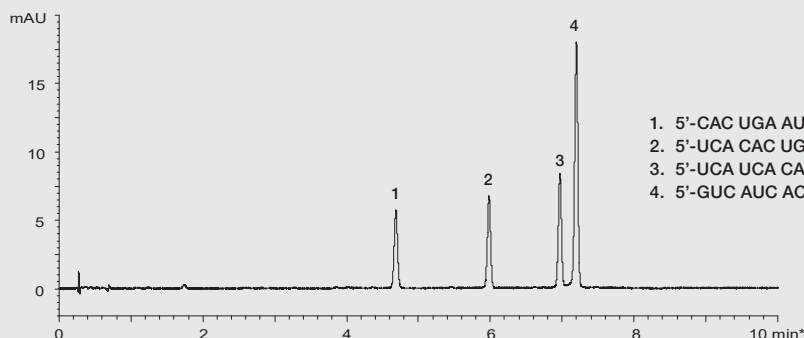


RP

Column: YMC-Triart C18 (1.9 μ m, 12 nm)
 50 x 2.1 mm ID
Part No.: TA12SP9-05Q1WT
Eluent: A) 200 mM HFIP*·8 mM triethylamine
 B) methanol

Gradient: 10–20%B (0–10 min)
Flow rate: 0.42 mL/min
Temperature: 65°C
Detection: UV at 260 nm
Injection: 1 μ L (2–4 nmol/mL)

LC/MS analysis of miRNA

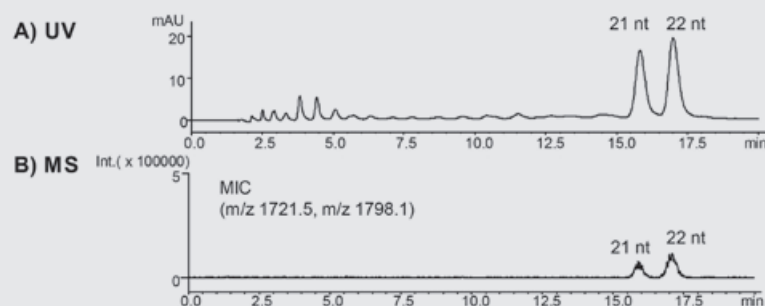

Courtesy of M. Yamada, SHIMADZU CORPORATION

5'-pUGG AGU GUG ACA AUG GUG UUG-3'
 5'-pUGG AGU GUG ACA AUG GUG UUG U-3'

(21mer, MW 6,890.1)
 (22mer, MW 7,196.3)


RP

Column: YMC-Triart C18 (3 μ m, 12 nm) 150 x 2.0 mm ID
Part No.: TA12S03-1502WT
Eluent: A) 10 mM di-n-butylamine-acetic acid (pH 7.5)
 B) 10 mM di-n-butylamine-acetic acid
 (pH 7.5)/acetonitrile (50/50)

Gradient: 62–72%B (0–20 min)
Flow rate: 0.2 mL/min
Temperature: 30°C
Detection: A) UV at 260 nm
 B) ESI-negative mode

Injection: 4 μ L (5 nmol/mL)
System: LC) Shimadzu Prominence
 MS) Shimadzu LCMS2020

Analysis of phosphoramidites

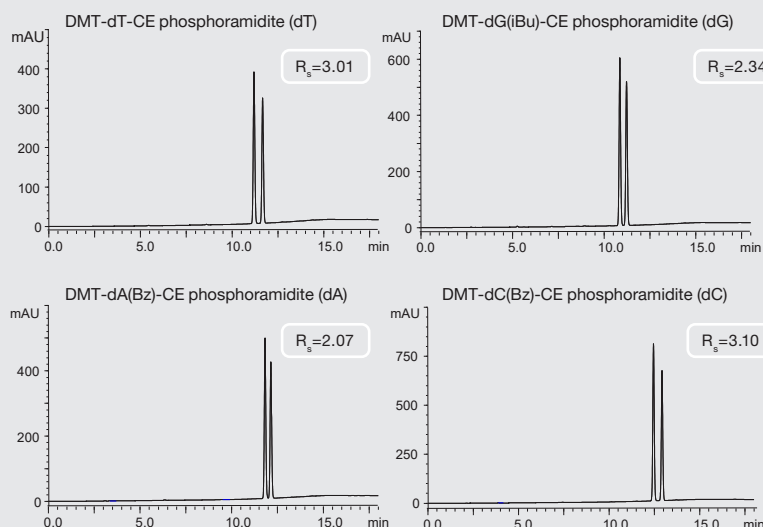

RP

Column: YMC-Triart C18
 (5 μ m, 12 nm)
 150 x 4.6 mm ID

Part No.: TA12S05-1546WT

Eluent: A) 20 mM triethylammonium acetate (pH 7.0) /
 acetonitrile (90/10)
 B) acetonitrile

Gradient: 45–95%B (0–12 min),
 95%B (12–17 min)

Flow rate: 1.0 mL/min

Temp.: 30°C

Detection: UV at 254 nm


Injection: 4 μ L (1.0 mg/mL)

OLIGONUCLEOTIDE APPLICATIONS


LC-HRMS analysis of the antisense oligonucleotide Mipomersen (Kynamro®)

RP


Total ion chromatogram

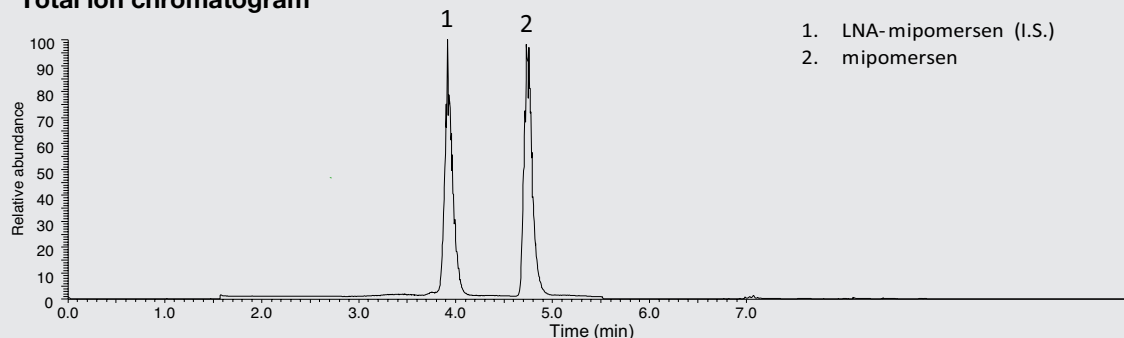

Mass spectrum of mipomersen

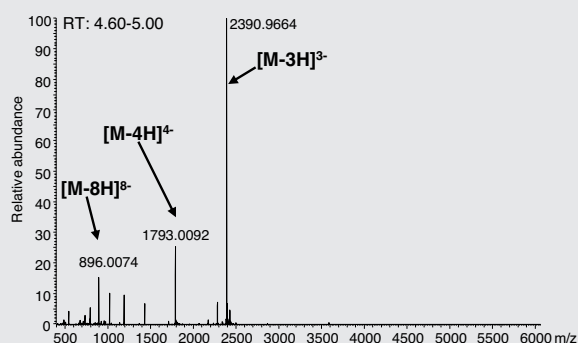

Mass spectrum of LNA-mipomersen

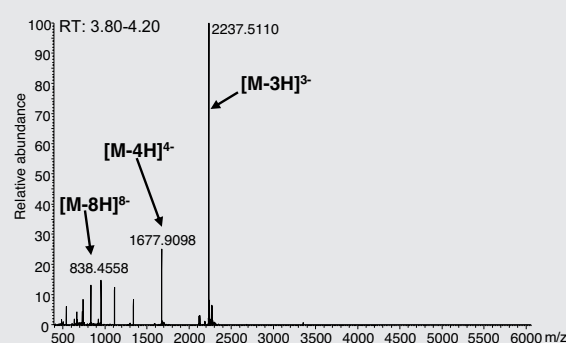

Column: YMC-Triart C8 metal-free PEEK-lined (1.9 μ m, 12 nm)[†] 100 x 2.1 mm ID
Part No.: T012SP9-10Q1PTP
Eluent: A) water/triethylamine/HFIP^{*2} (100/0.4/2; triethylamine 28.0 mM, HFIP 135.8 mM)
 B) methanol/triethylamine/HFIP (100/0.4/2)
Gradient: [Sample separation step]
 10–40%B (0–5.0 min)
 [Column wash steps]
 40–70%B (5.0–5.1 min), 70%B (5.1–7.0 min), 70–10%B (7.0–7.1 min), 10%B (7.1–8.0 min), 10–90%B (8.0–8.1 min), 90%B (8.1–9.0 min),
 90–10%B (9.0–9.1 min), 10%B (9.1–10.0 min), 10–90%B (10.0–10.1 min), 90%B (10.1–11.0 min), 90–10%B (11.0–11.1 min)
Flow rate: 0.3 mL/min
Temperature: 50°C
Injection: 10 μ L (1000 ng/mL)
System: LC) Vanquish Binary Pump H system
 HRMS) Orbitrap HRMS Q Exactive Plus

*1 Prewash the column prior to the first use with water/methanol/phosphoric acid (70/30/0.1) for 1 hour


*2 1,1,1,3,3,3-hexafluoro-2-propanol

Reference: Y. Sun et al, Development of a bioanalytical method for an antisense therapeutic using high-resolution mass spectrometry, Bioanalysis, 2020 NOV 26, doi: 10.4155/bio-2020-0225.

OLIGONUCLEOTIDE APPLICATIONS

mRNA coupled to enhanced green fluorescent protein (EGFP)

RP

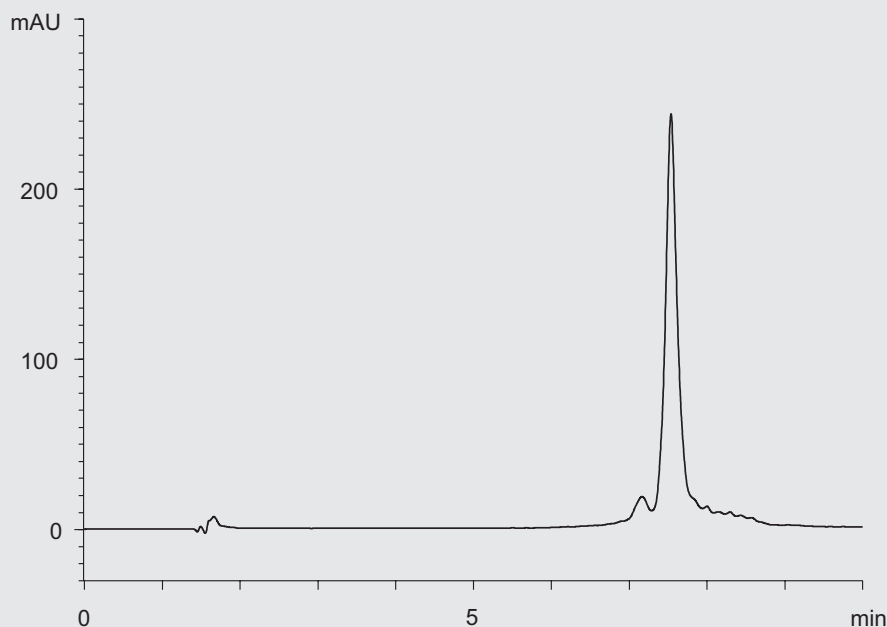

Column: YMC Accura Triart Bio C4
(3 μ m, 30 nm)
100 x 2.1 mm ID

Part No.: TB30S03-10Q1PTC

Eluent: A) 50 mM TEAA* (pH 7.0)
acetonitrile (95/5)
B) acetonitrile

Gradient: 5–10%B (0–10 min)

Flow rate: 0.2 mL/min

Temperature: 80 °C

Detection: UV at 254 nm


Injection: 2 μ L (0.25 mg/mL)

Sample: CleanCap® EGFP mRNA
(5 moU) (TriLink Bio
Technologies)

*triethylammonium acetate

IEX analysis of EGFP mRNA (996mer)

AEX

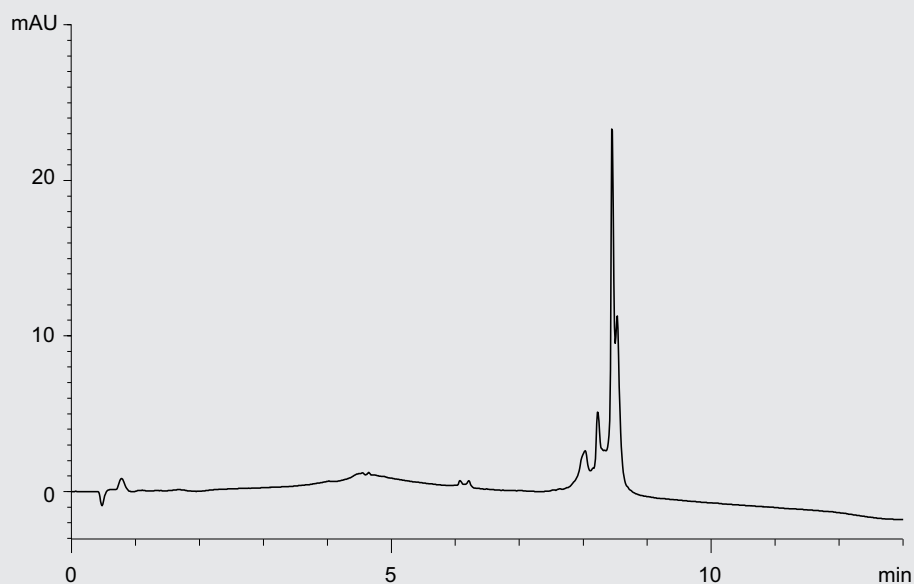

Column: BioPro IEX QF (5 μ m)
100 x 4.6 mm ID

Part No.: QF00S05-1046WP

Eluent: A) 10 mM NaOH (pH 12)
B) 10 mM NaOH (pH 12)
containing 2M NaCl

Gradient: 0–100%B (0–9 min),
100% B (9–13 min)

Flow rate: 1.0 mL/min

Temperature: 15 °C

Detection: UV at 260 nm


Injection: 5 μ L (0.025 mg/mL)

Sample: CleanCap® EGFP mRNA
(5 moU) (TriLink Bio
Technologies)

OLIGONUCLEOTIDE APPLICATIONS

siRNA duplex under denaturing conditions

RP

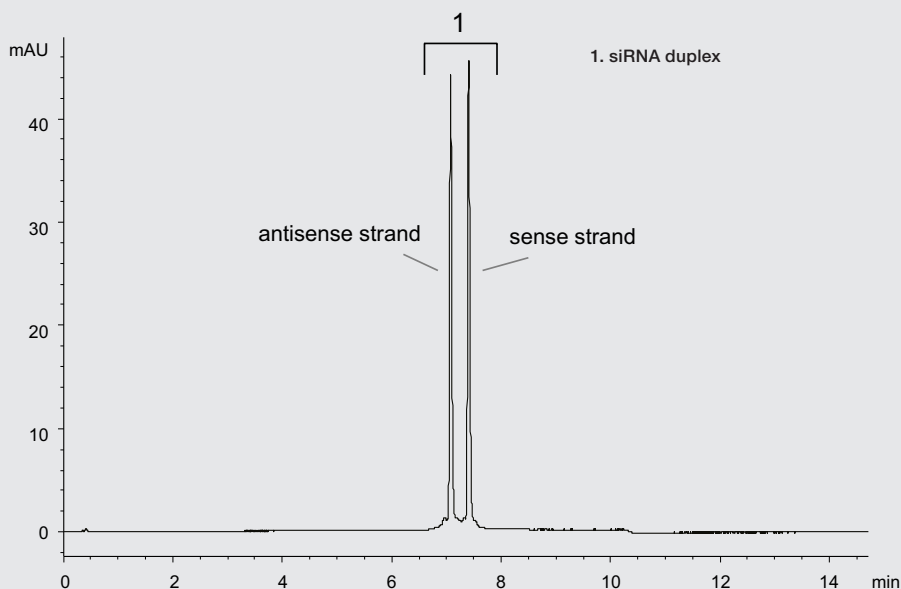

Column: YMC Accura Triart Bio C18 (1.9 μ m, 30 nm) 50 x 2.1 mm ID

Part No.: TA30SP9-05Q1PTC

Eluent: A) 15 mM TEAA* (pH 8)
B) methanol

Gradient: 5%–20%B (0–15 min)

Flow rate: 0.42 mL/min

Temperature: 65 °C

Detection: UV at 260 nm


Injection: 1 μ L (5 nmol/mL)

Sample: siRNA duplex

*triethylammonium acetate

siRNA duplex under denaturing conditions

AEX

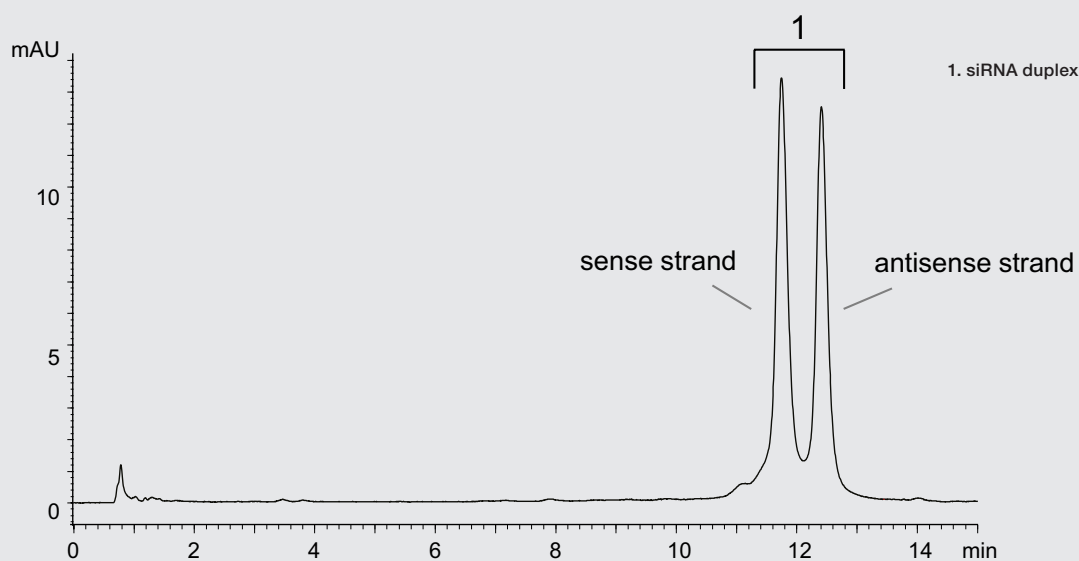

Column: BioPro IEX QF (5 μ m) 100 x 4.6 mm ID

Part No.: QF00S05-1046WP

Eluent: A) 10 mM NaOH
B) 10 mM NaOH containing 1 M NaClO₄

Gradient: 30%–37%B (0–15 min)

Flow rate: 1.0 mL/min

Temperature: 25 °C


Detection: UV at 260 nm

Injection: 4 μ L (5 nmol/mL)

Sample: siRNA duplex

OLIGONUCLEOTIDE APPLICATIONS

Challenging phosphorothioate oligonucleotides

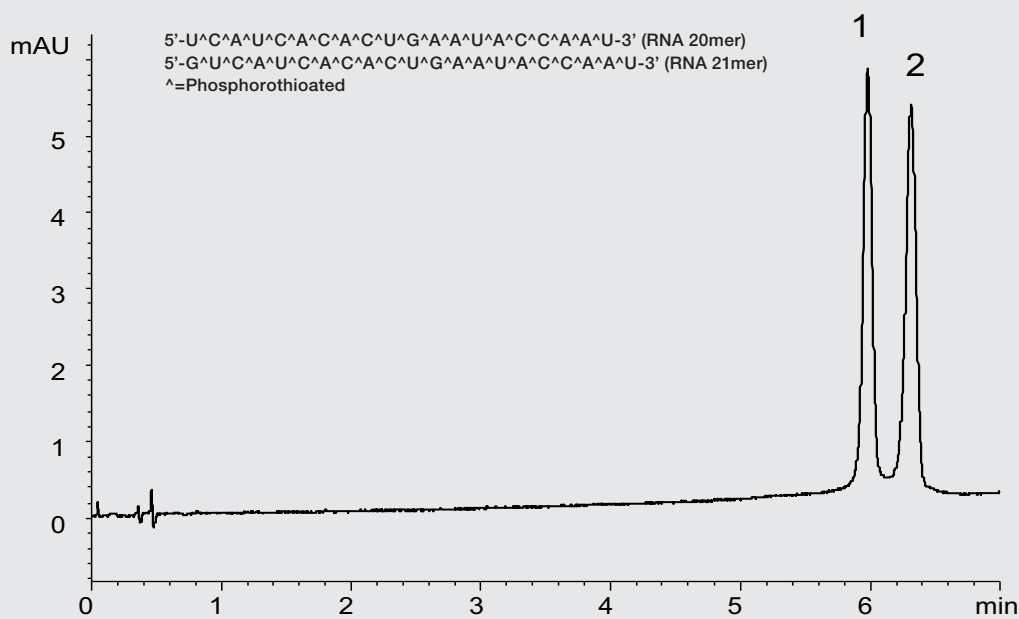


RP

Column: YMC Accura Triart Bio C18
(1.9 µm, 30 nm) 50 x 2.1 mm ID
Part No.: TA30SP9-05Q1PTC
Eluent: A) 15 mM triethylamine - 400 mM HFIP*
B) methanol

Gradient: 10–20%B (0–10 min)
Flow rate: 0.42 mL/min
Temperature: 65°C
Detection: UV at 260 nm
Injection: 1 µL (each 1.0 nmol/mL)

* 1,1,1,3,3,3-hexafluoro-2-propanol

AEX analysis of siRNA duplex under non-denaturing conditions

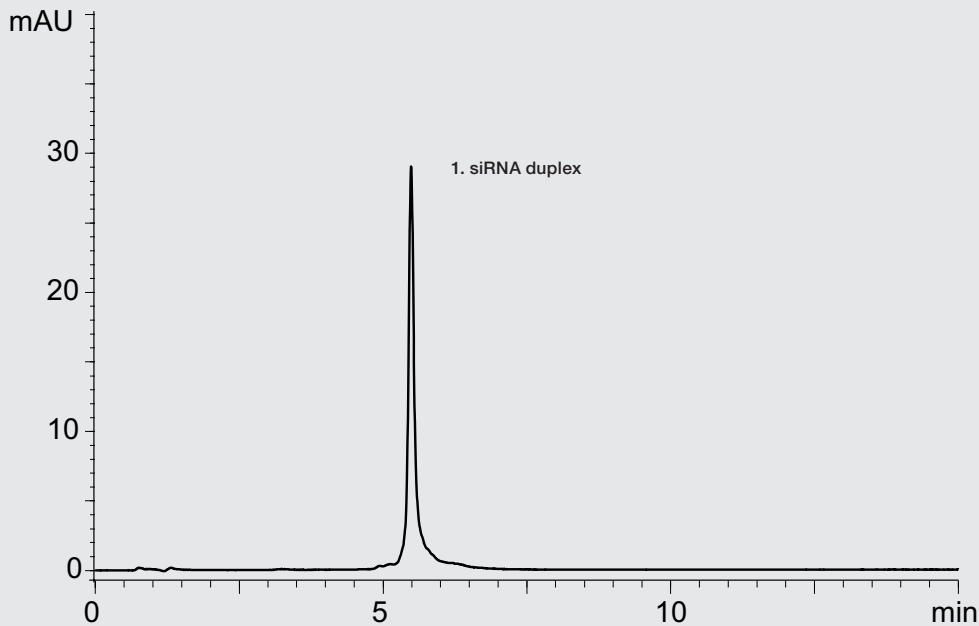

AEX



Column: BioPro IEX QF (5 µm) 100 x 4.6 mm ID
Part No.: QF00S05-1046WP
Eluent: A) 20 mM Tris-HCl (pH 8.1)
B) 20 mM Tris-HCl (pH 8.1) containing 1 M NaClO₄

Gradient: 25%–40%B (0–15 min)
Flow rate: 1.0 mL/min
Temperature: 25°C
Detection: UV at 260 nm
Injection: 4 µL (5 nmol/mL)

Sample: siRNA duplex

YMC IP-RP

YMC COLUMN

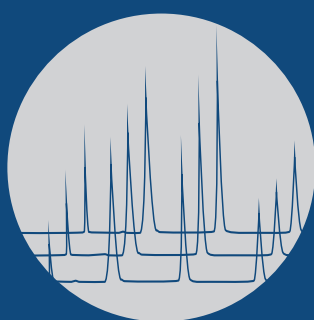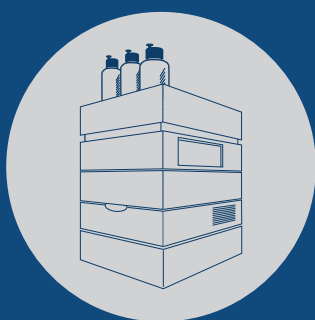



IP-RP – UHPLC/HPLC SELECTIVITIES

FEATURES

- Selection of C18, C8 and C4 columns
- For ion-pairing (IP-RP) UHPLC and HPLC
- pH- and temperature stable phases
- Superior reproducibility
- Bioinert coated YMC Accura column hardware

Selectivities for oligonucleotides

	YMC-Triart C18	YMC-Triart Bio C18	YMC-Triart C8	YMC-Triart Bio C4	Hydrosphere C18
Base particle	organic/inorganic hybrid silica				silica
Modification	C18 (USP L1)	C18 (USP L1)	C8 (USP L7)	C4 (USP L26)	C18 (USP L1)
Particle size / μm	1.9, 3, 5	1.9, 3, 5	1.9, 3, 5	1.9, 3, 5	2, 3, 5
Pore size / nm	12	30	12	30	12
pH range	1.0 – 12.0	1.0–12.0	1.0–12.0	1.0–10.0	2.0–8.0
Temp. Range	pH < 7: 90 °C pH > 7: 50 °C	pH < 9: 90 °C pH > 9: 50 °C	pH < 7: 90 °C pH > 7: 50 °C	pH < 7: 90 °C pH > 7: 50 °C	50 °C
Column hardware	bioinert coated YMC Accura				stainless-steel

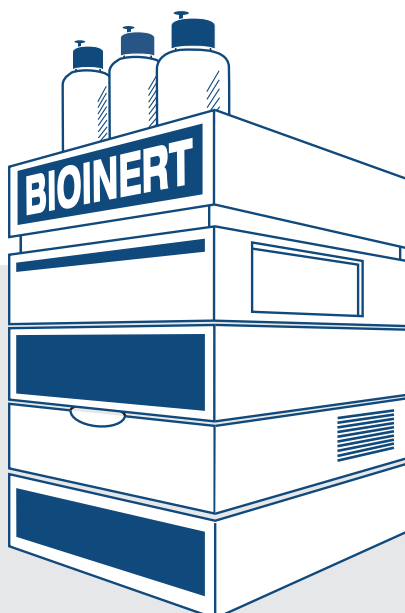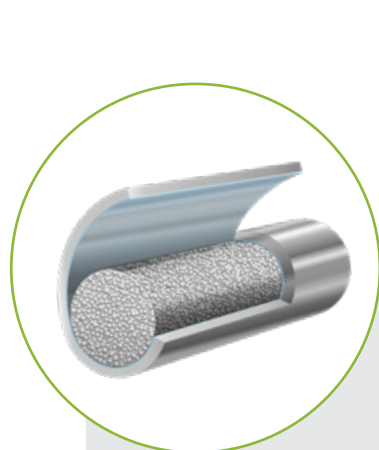


**Bioinert
hardware
available!**

Bioinert YMC Accura Triart columns are available for improved sensitivity and peak shape of coordinating compounds such as nucleotides or oligonucleotides.


IP-RP – UHPLC/HPLC SELECTIVITIES

Column Selection Tool according to nucleotide length

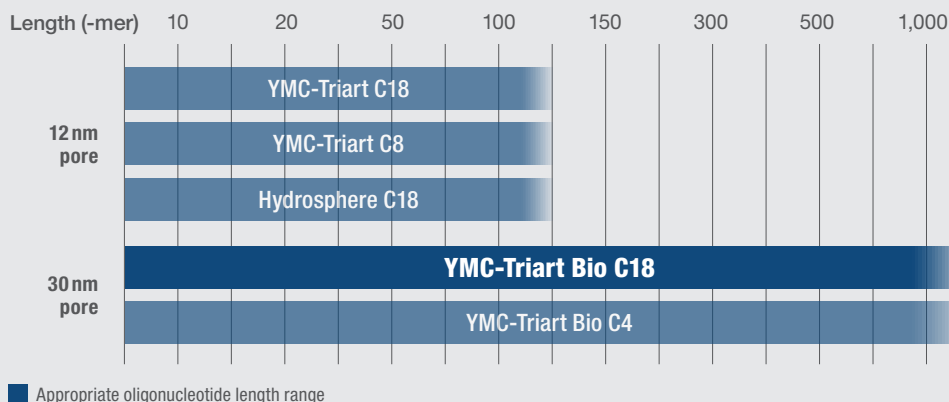

First choice for oligonucleotides of various lengths

Ideal choice for challenging analytes such as phosphorothioate oligonucleotides

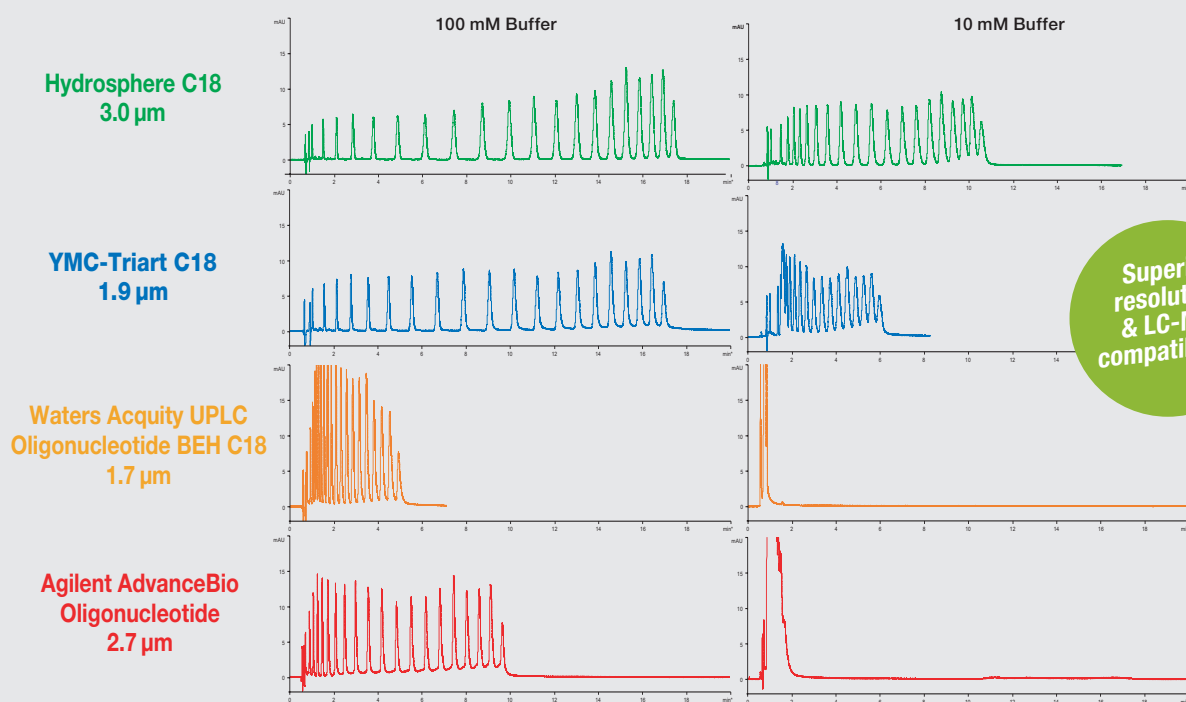


Column: 50 x 2.1 mm ID
Eluent: A) 100 mM or 10 mM triethylamine-acetic acid (pH 6)
 B) A/acetonitrile (80/20)
Gradient: 50–65%B (0–20 min)

Flow rate: 0.21 mL/min
Temperature: 35 °C
Detection: UV at 260 nm
Injection: 2 µL (5 nmol/mL)
Sample: Oligodeoxythymidylic acid [d(pT)2-20]

The separation of oligo(deoxythymidylic acids), d(pT)2-20, was compared using 100 mM or 10 mM triethylammonium acetate (TEAA) buffer, under the same gradient conditions. Both Hydrosphere C18 and YMC-Triart C18 showed enhanced retention and resolution compared to other commercially available C18 phases designed for oligonucleotide analysis, even at the low ion-pairing buffer concentration such as 10 mM. The higher concentration provides stronger retention and superior resolution of oligonucleotides, although a lower concentration has the advantages of increasing the signal intensity and reducing system contamination in HPLC-MS analysis.

IP-RP – UHPLC/HPLC SELECTIVITIES

Optimum resolution for short and long poly(dT) oligonucleotides

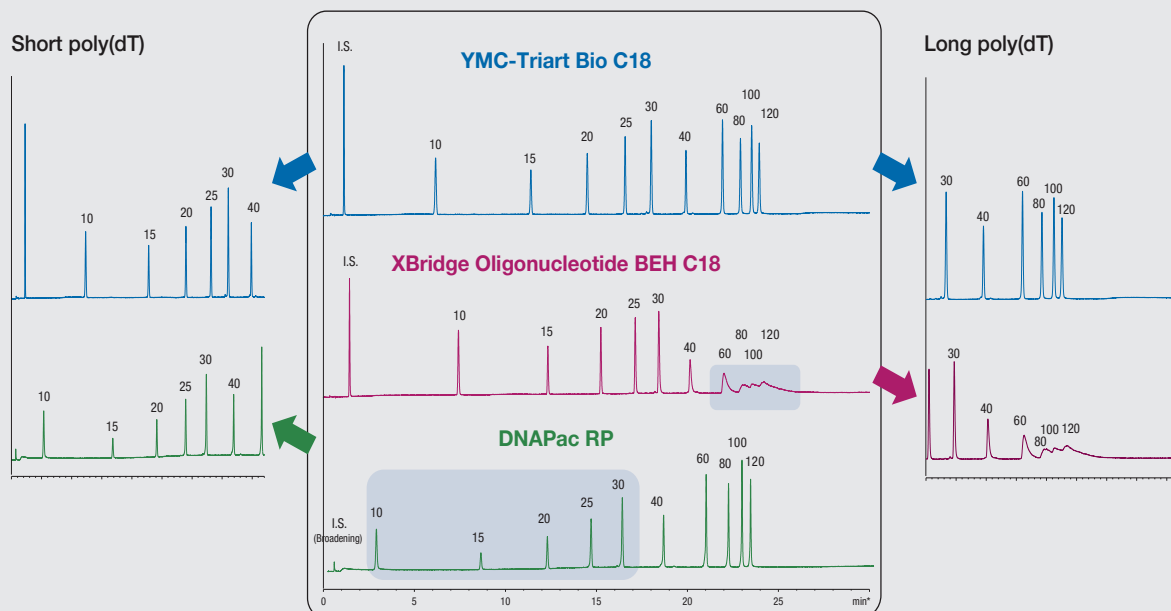

Columns: YMC-Triart Bio C18 (3 μ m, 30 nm)
50 x 2.1 mm ID
XBridge Oligonucleotide BEH C18
(2.5 μ m, 13 nm) 50 x 2.1 mm ID
DNAPac RP (4 μ m, proprietary)
50 x 2.1 mm ID

Part No.: TA30S03-05Q1WT
Eluent: A) 4 mM TEA* - 100 mM HFIP**
B) methanol
Gradient: 0.5%B/min, initial %B=5
Flow rate: 0.42 mL/min
Detection: UV at 260 nm


Temperature: 65 °C
Injection: 1.0 μ L
Sample: Poly(dT) oligonucleotides

*triethylamine, **1,1,1,3,3,3-hexafluoro-2-propanol

The YMC-Triart Bio C18 column demonstrates a better resolution, higher recovery and reproducibility of poly(dT) oligonucleotides compared to the other two competitor columns. Longer poly(dT) oligonucleotides (60–120mer) were separated poorly by the competitor's hybrid silica based column, whereas YMC-Triart showed high resolution for oligonucleotides of all sizes. Peak areas and therefore recoveries of shorter poly(dT) oligonucleotides (10–40mer) were much smaller when separated using the dedicated DNA competitor column. In addition, YMC-Triart Bio C18 showed reproducible behaviour such as consistent peak areas. This makes YMC-Triart Bio C18 an ideal tool for analysis of poly(dT) oligonucleotides.


IP-RP – TEMP. & MOBILE PHASE EFFECTS

Phosphate buffer (pH 11.5, 40 °C)

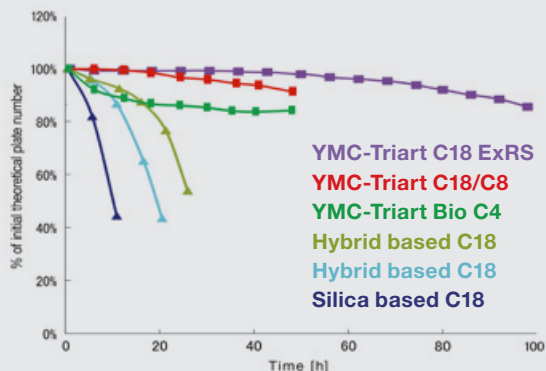


Column: 5 μ m, 150 x 4.6 mm ID
Eluent: 50 mM K_2HPO_4 - K_3PO_4 (pH 11.5)/methanol (90/10)
Flow rate: 1.0 mL/min
Temperature: 40 °C
Sample: benzyl alcohol

15mM TEA - 400mM HFIP (pH 8, 90 °C)

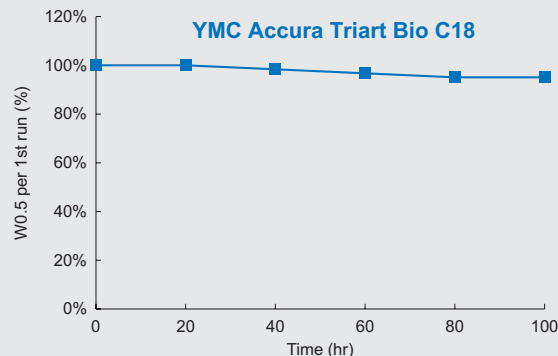

Column: YMC Accura Triart Bio C18 (1.9 μ m, 30 nm), 50 x 2.1 mm ID
Eluent: A) 15 mM TEA-400 mM HFIP (pH 8)
 B) methanol
Gradient: 6–16%B (0–10 min)
Flowrate: 0.42 mL/min
Temperature: 90 °C
Sample: 20mer RNA

Effect of four different ion pair reagents on the denaturation of a siRNA duplex at different temperatures

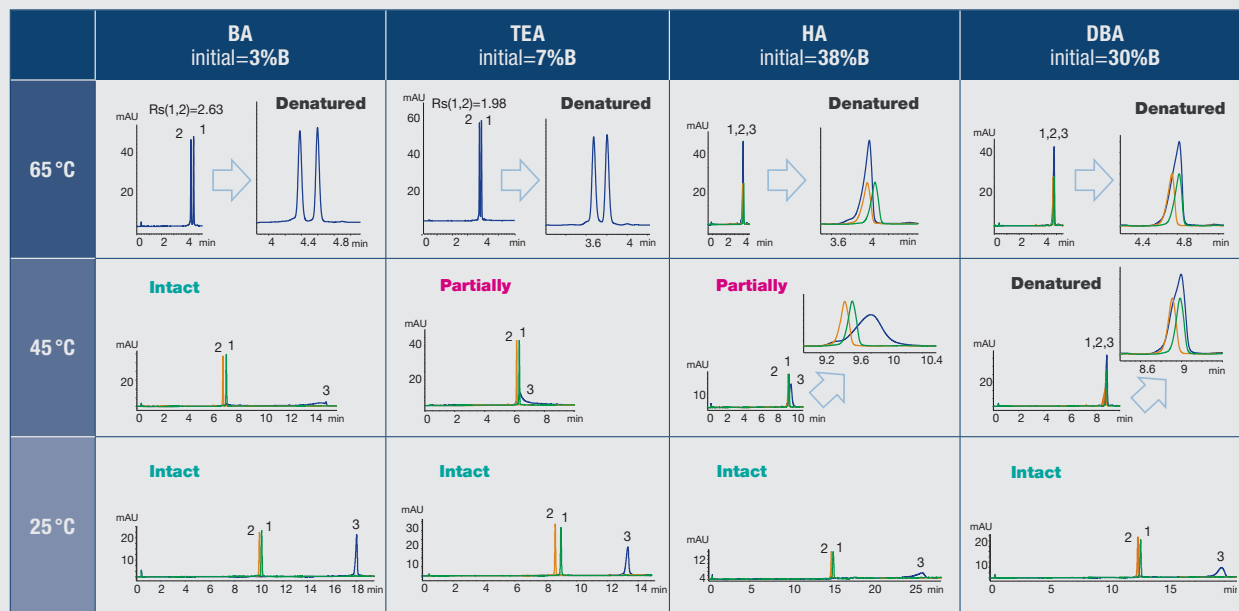

1. sense strand 2. antisense strand 3. siRNA duplex

Column: YMC Accura Triart Bio C18 (1.9 μ m, 30 nm), 50 x 2.1 mm ID
Part No.: TA30SP9-05Q1PTC
Eluent: A) 8 mM alkylamine - 100 mM HFIP* (pH 8)
 B) methanol
Gradient: 1%B/min
Flow rate: 0.42 mL/min


Temperature: 65 °C, 45 °C or 25 °C
Detection: UV at 260 nm
Injection: 1 μ L (5 nmol/mL)
Sample: siRNA duplex, sense strand, antisense strand
 5'-CGU ACG CGG AAU ACU UCG AdTdT-3' sense strand
 3'-dTdTGCA UGC GCC UUA UGA AGC U-5' antisense strand

*1,1,1,3,3,3-hexafluoro-2-propanol

IP-RP – EXPERT TIPS: ION-PAIRING SALTS

Effect of composition and salt concentration of ion-pairing mobile phase on the separation and signal intensity

Comparison of separation and ESI-MS signal intensity using different ion-pairing buffers and organic solvents

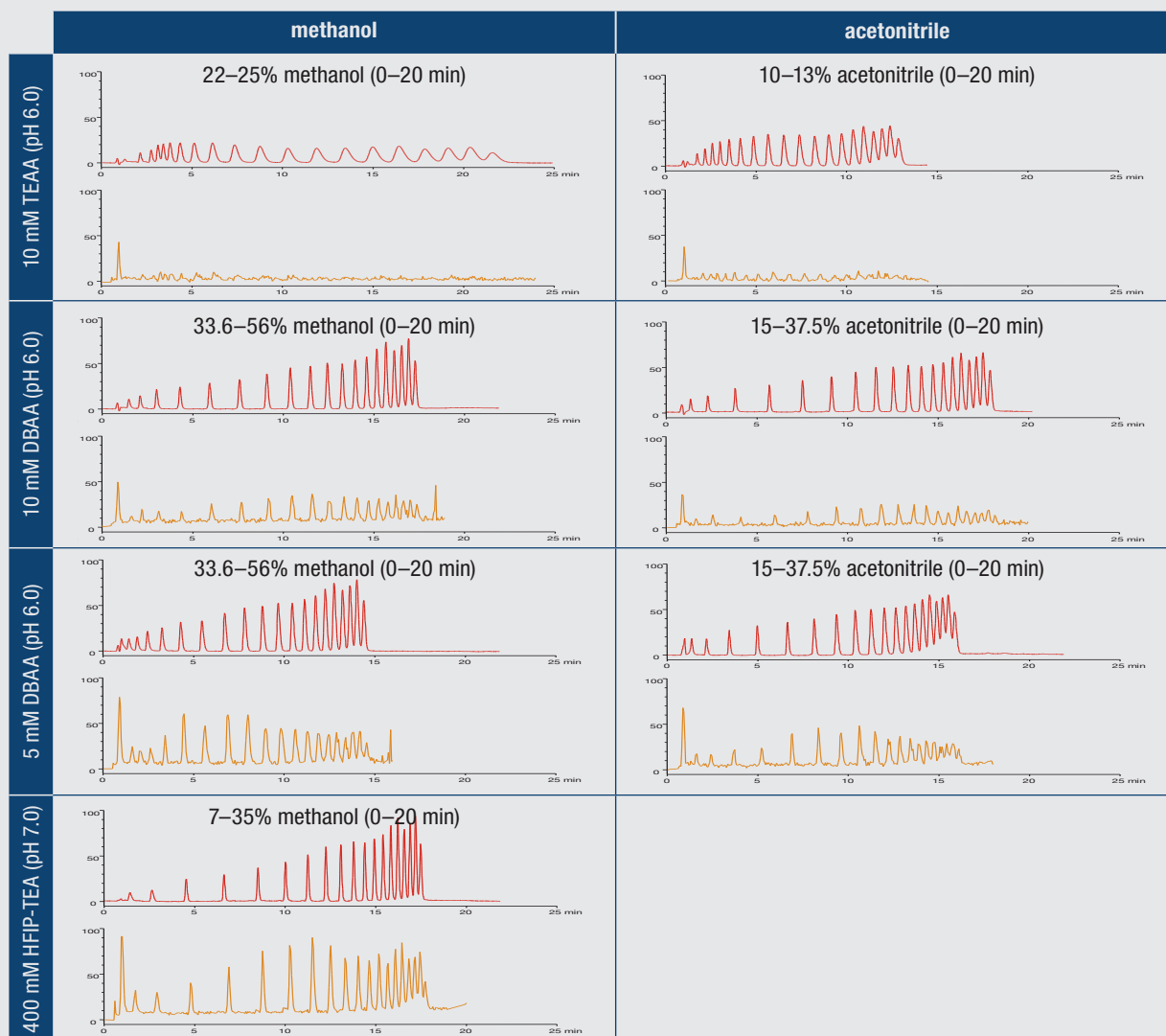

Columns: Hydrosphere C18 (3.0 μ m, 12 nm)
50 x 2.0 mm ID
Part No.: HS12S03-0502WT
Flow rate: 0.2 mL/min
Temperature: 35°C

Detection: UV at 269 nm (red), ESI negative mode (orange)
Injection: 5 μ L (25 pmol/mL)
Sample: Oligo(deoxythymidylic acid) [dp (dT) 2-20]


eluent abbreviations
TEAA: triethylammonium acetate
DBAA: di-n-butylammonium acetate
HFIP: 1,1,1,3,3,3-hexafluoro-2-propanol
TEA: triethylamine

The mobile phase composition has different effects on the separation and signal intensity in electrospray ionisation mass spectrometry (ESI-MS) of oligonucleotides. Using different gradient conditions, acceptable retention and resolution can be achieved (upper UV chromatograms; red trace) for each separation by optimising the gradient slope of the organic solvent regardless of the type of mobile phase. The ESI-MS intensity is significantly influenced by the type and concentration of ion-pairing

buffer as shown in the lower MS chromatograms (orange trace). HFIP-TEA buffer/methanol systems provide the maximum MS intensity. Enhanced retention and MS intensity are obtained using 10 mM DBAA buffer compared to 10 mM TEAA buffer, and the lower DBAA concentration results in approximately 1.5–3 times increase in the intensity without any change in the concentration of organic solvent.

YMC ACCURA BIOINERT HARDWARE


YMC COLUMN

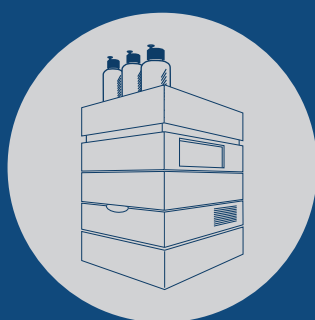

YMC ACCURA BIOINERT HARDWARE


BIOINERT COLUMNS FOR BIOSEPARATIONS AND COORDINATING COMPOUNDS

- Exceptional peak shapes with high sensitivities
- Excellent recoveries without column preconditioning
- Superior reproducibility and no carry-over effects
- Ideal for highly sensitive LC/MS analyses

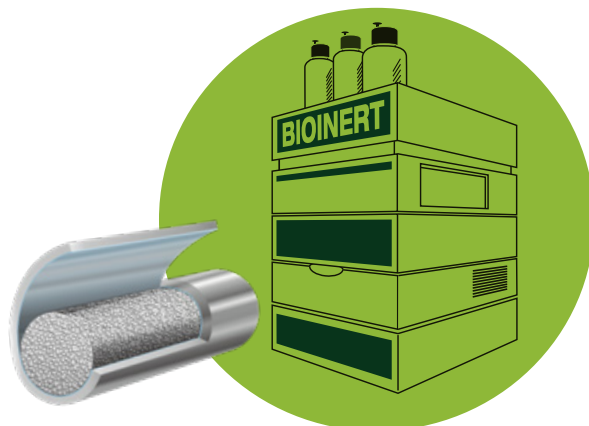

Specifications


	YMC Accura Triart	YMC Accura BioPro IEX
Modifications	C18, Bio C18, C8, Bio C4, Diol-HILIC	QF
Particle Size	1.9, 3 and 5 μm	3 and 5 μm
Column hardware	stainless-steel with bioinert coating	
Frit hardware	stainless-steel with bioinert coating	
Pressure limit	1.9 μm : 100 MPa (15,000 psi) 3/5 μm : 45 MPa (6,525 psi)	3 μm : 15–30 MPa (2,175–4,350 psi) 5 μm : 10–30 MPa (1,450–4,350 psi)

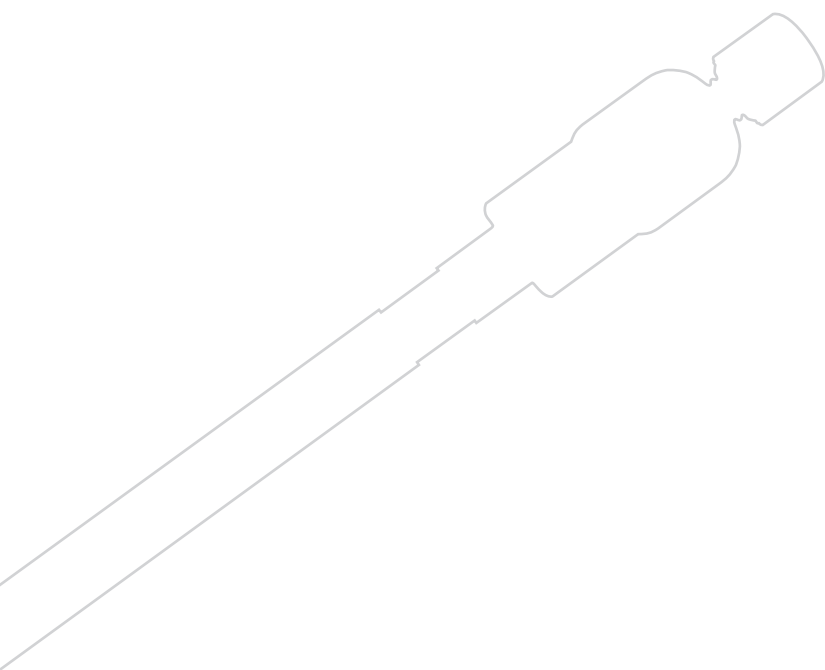


YMC ACCURA BIOINERT HARDWARE

Improved sensitivity for coordination compounds


1. ATP

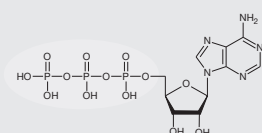


2. ADP

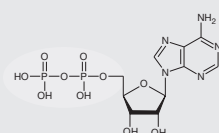


3. AMP

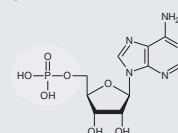

YMC Accura Triart C18

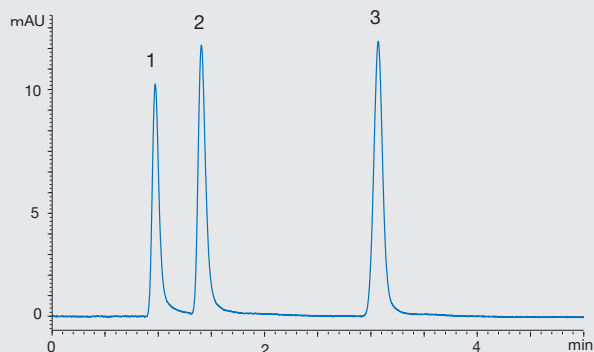

Standard column

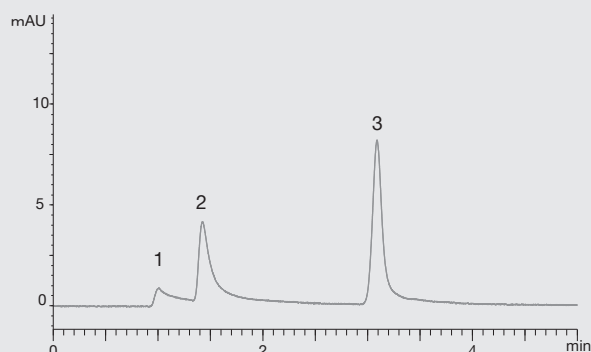


Column: YMC Accura Triart C18 (1.9 μ m, 12 nm) 50 x 2.1 mm ID (bioinert hardware)
YMC-Triart C18 (1.9 μ m, 12 nm) 50 x 2.1 mm ID (standard hardware)
Part Nos.: TA12SP9-05Q1PTC
TA12SP9-05Q1WT
Eluent: 5 mM HCOONH₄

Flow rate: 0.21 mL/min
Temperature: 25 °C
Detection: UV at 265 nm
Injection: 1 μ L (10 μ g/mL)
System: bioinert/"metal-free" HPLC system

Metal coordinating compounds, which have a phosphate group in their structure, tend to show poor peak shape due to interactions with metals, such as the stainless-steel in column bodies and frits. By using a bioinert column hardware, better peak shapes can be expected. Nucleotides with phosphate groups also show better peak shapes when compared to the regular column hardware. The bioinert coated YMC Accura column hardware is ideal for highly sensitive analyses using LC/MS.

YMC ACCURA BIOINERT HARDWARE


YMC ACCURA: DURABLE BIOINERT COATING

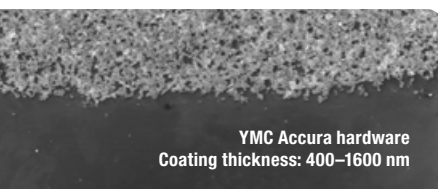

The robust bioinert coating used on YMC Accura hardware is 130 to 320-fold thicker, making it more durable than other similar hardware concepts. A long-term inertness against sensitive substances is ensured.


In order to demonstrate its robustness, a YMC Accura column was packed multiple times. Even though this process is hard on the column surface, the coating remains unaffected (SEM* picture: top area is bare steel for comparison).

*Scanning Electron Microscope

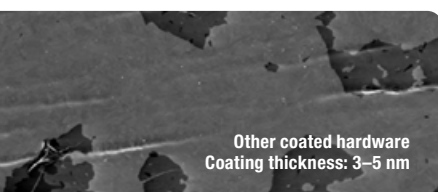

Other coated columns can lose their inertness over time. This will again lead to adsorption of sensitive compounds on the uncovered metallic surfaces. Peak tailing, loss of recovery and sample carry-over are typical results of the delamination of the coating. On a competitor's column, most of the coating was already delaminated after merely unpacking the column. (The dark spots in this picture are the remaining coating.)

High surface inertness without any adsorption

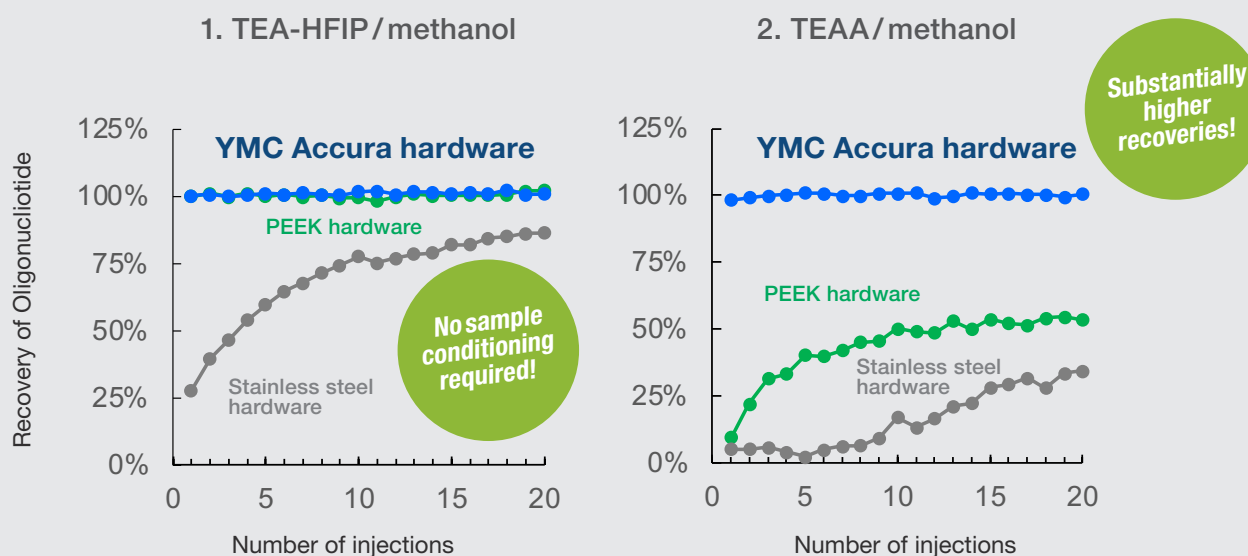

Column: Empty YMC Accura (50 x 2.1 mm ID, w/o stationary phase)
Eluent: 1) 8 mM TEA 200 mM HFIP / methanol (82/18)
 2) 100 mM TEAA / methanol (82/18)
Flow rate: 0.42 mL/min
Detection: UV at 260 nm

Temperature: 65°C
Injection: 1 µL (18 nmol/mL)
Sample: All PS RNA 20mer (1) (5'-U[△]C[△]A[△]U[△]C[△]A[△]C[△]A[△]C[△]U[△]G[△]A[△]A[△]U[△]A[△]C[△]C[△]A[△]A[△]U-3')


[△]=Phosphorothioate

The YMC Accura hardware with its inert surface area prevents adsorption of oligonucleotides using a range of different buffers. No sample conditioning is required. YMC Accura columns further provide significantly higher recoveries and sensitivities that cannot be achieved with regular stainless-steel columns – even after conditioning with 20 sample injections. These ready-to-use columns ensure high recovery and reproducibility from the very first use.

YMC ACCURA BIOINERT HARDWARE

SIGNIFICANTLY HIGHER SENSITIVITY AND RECOVERY

Ideal choice for challenging analytes such as phosphorothioate oligonucleotides

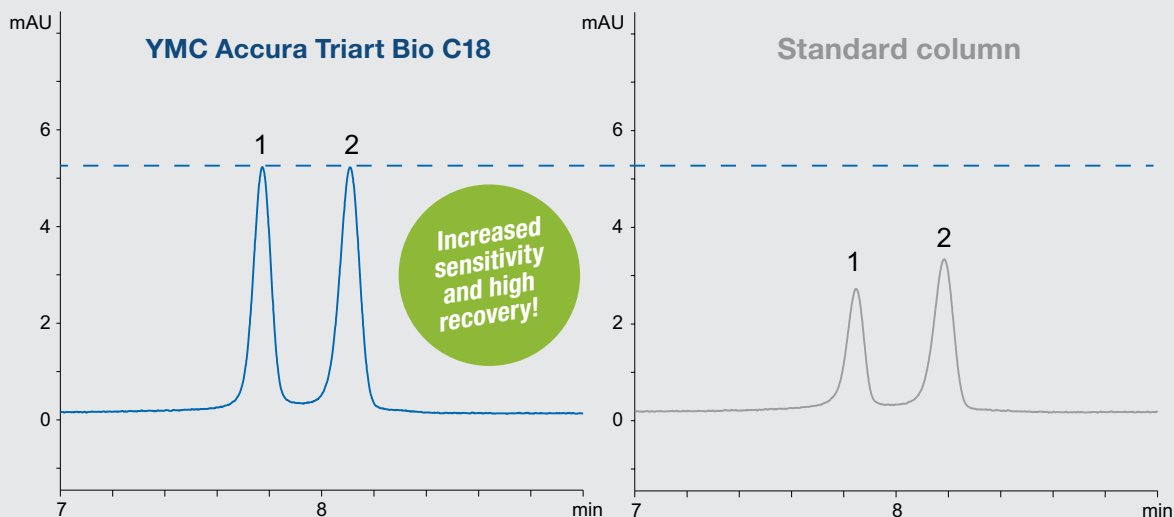


Column: YMC Accura Triart Bio C18 (1.9 µm, 30 nm) 50 x 2.1 mm ID
Part No.: TA30SP9-05Q1PTC
Eluent: A) 15 mM triethylamine - 400 mM HFIP*
 B) methanol
Gradient: 8–18%B (0–10 min)
Flow rate: 0.42 mL/min
Temperature: 65°C

Detection: UV at 260 nm
Injection: 1 µL
Sample: All PS RNA 20mer (1) (5'-U[^]C[^]A[^]A[^]U[^]C[^]A[^]C[^]A[^]C[^]U[^]G[^]A[^]A[^]U[^]A[^]C[^]A[^]A[^]U[^]-3')
 All PS RNA 21mer (2) (5'-G[^]U[^]C[^]A[^]A[^]U[^]C[^]A[^]C[^]A[^]C[^]U[^]G[^]A[^]A[^]U[^]A[^]C[^]A[^]C[^]A[^]A[^]U[^]-3')
[^]=Phosphorothioate

*1,1,1,3,3,3-hexafluoro-2-propanol


High sensitivity and recovery

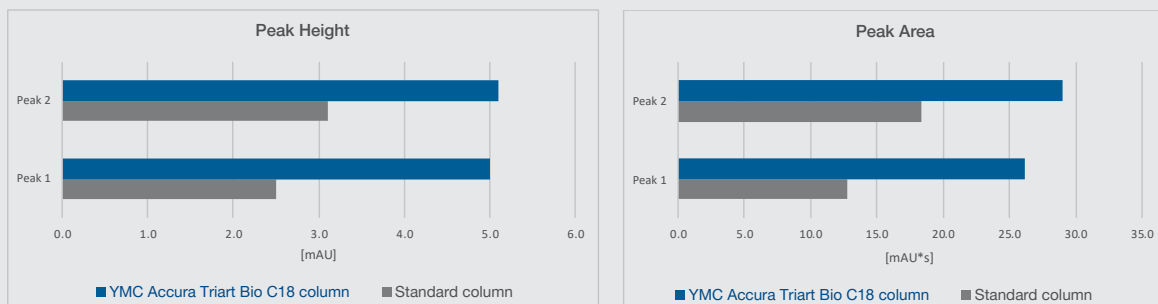


The YMC Accura Triart Bio C18 column provides double peak heights and peak areas for the oligonucleotides compared to those for regular stainless-steel columns. YMC Accura Triart columns enhance the analytical sensitivity significantly, thereby avoiding the loss of high-value samples.

YMC ACCURA BIOINERT HARDWARE

RELIABLE RESULTS FROM THE FIRST INJECTION

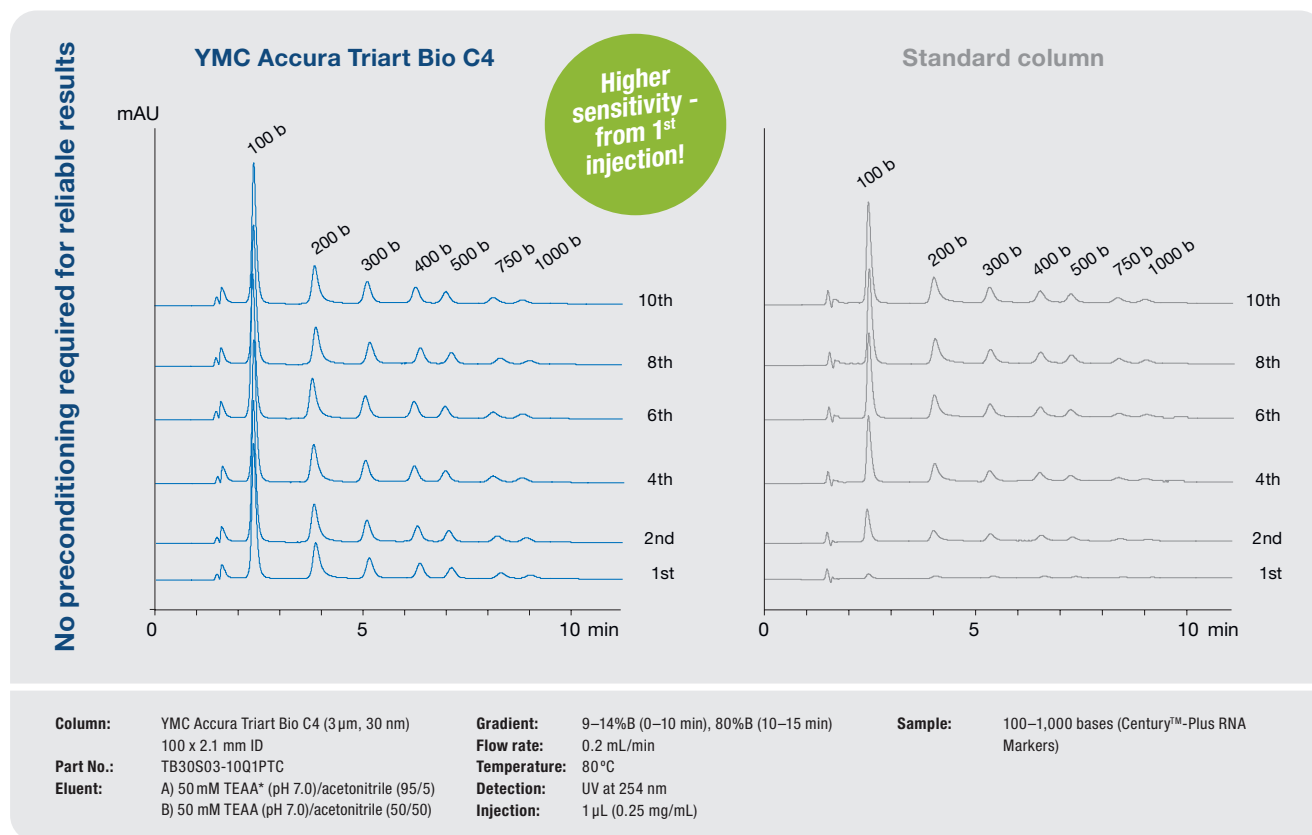



* Triethylammonium acetate

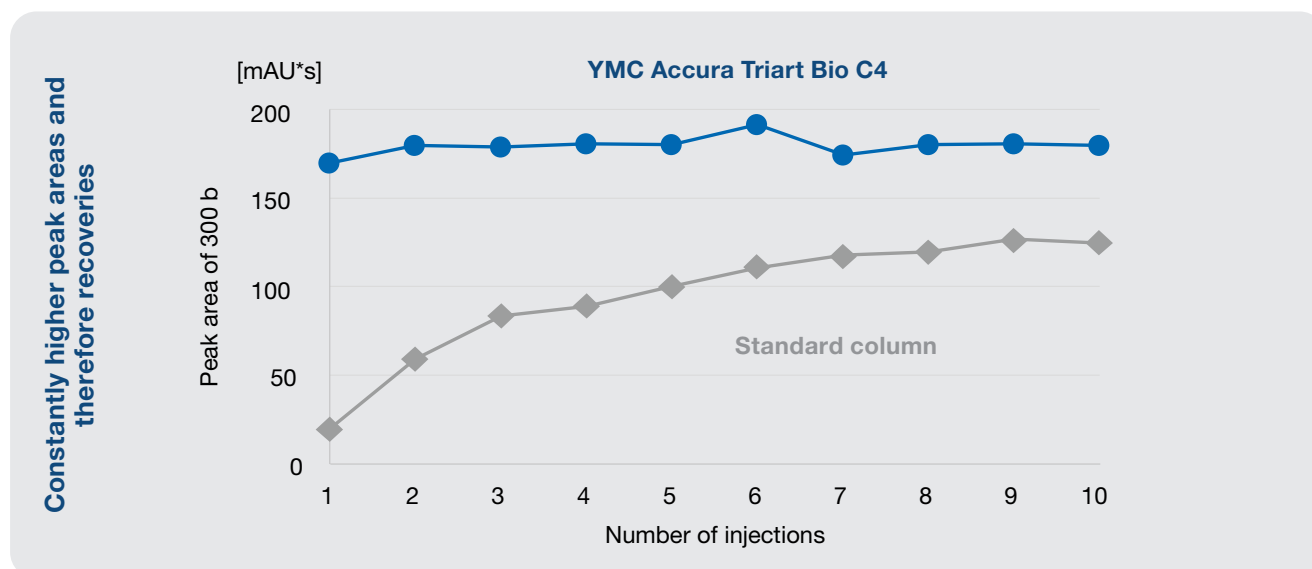

The YMC Accura Triart Bio C4 column shows stable peak areas from the first injection, while the standard stainless-steel column provides only 10% of the peak area (for the 300 base marker) with the first injection. Even after the tenth injection, the peak areas of the stainless-steel column are considerably less than those of the YMC Accura Triart column.

YMC AEX


YMC COLUMN

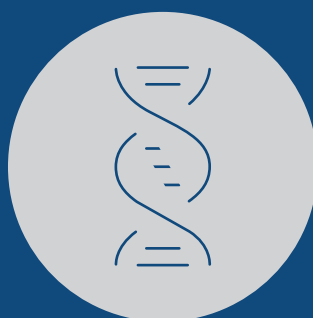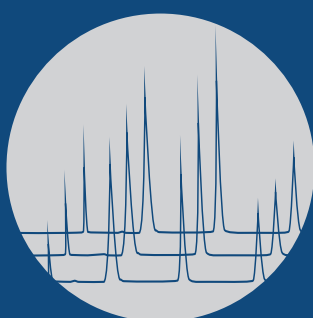


AEX – HPLC SELECTIVITY

FEATURES

- Non-porous hydrophilic polymers
- High recovery and high resolution
- Minimised nonspecific adsorption
- Excellent reproducibility
- Bioinert coated YMC Accura column hardware

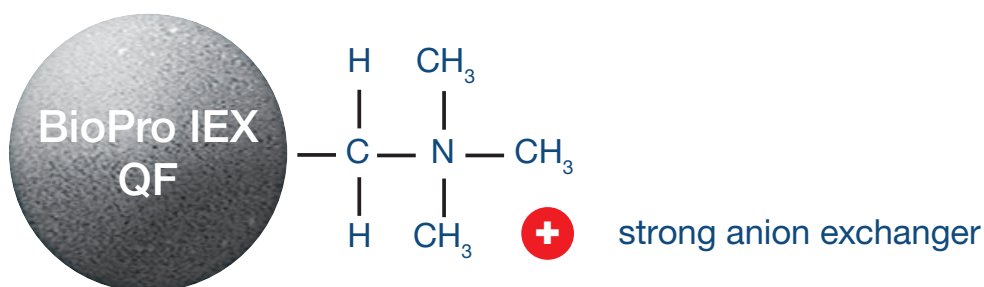


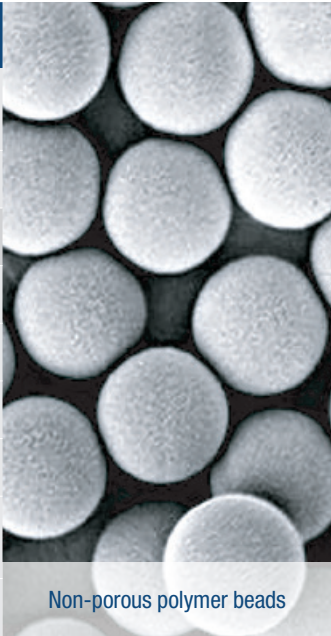	BioPro IEX QF	
Matrix	hydrophilic polymer (polymethacrylate)	
Particle size / μm	3, 5	
Pore size / nm	non-porous	
Charged group	$-\text{CH}_2\text{N}^+(\text{CH}_3)_3$	
Counter ion	Cl^-	
Available pH range	2.0–12.0	
Temperature range	4–80°C	
Pressure limit	3 μm : 15–30 MPa (2,175–4,350 psi) 5 μm : 10–30 MPa (1,450–4,350 psi)	
Column hardware	bioinert coated YMC Accura	Non-porous polymer beads

YMC's anion exchanger (AEX) columns of the BioPro IEX series are available with strong exchanger modification, based on 3 or 5 μm non-porous (QF columns) hydrophilic polymer beads. The non-porous particles offer high efficiency, very high resolution and low operating pressures.

AEX – LOADABILITY


SHARP PEAKS AND RELIABLE RECOVERY WITH DIFFERENT COLUMN DIMENSIONS

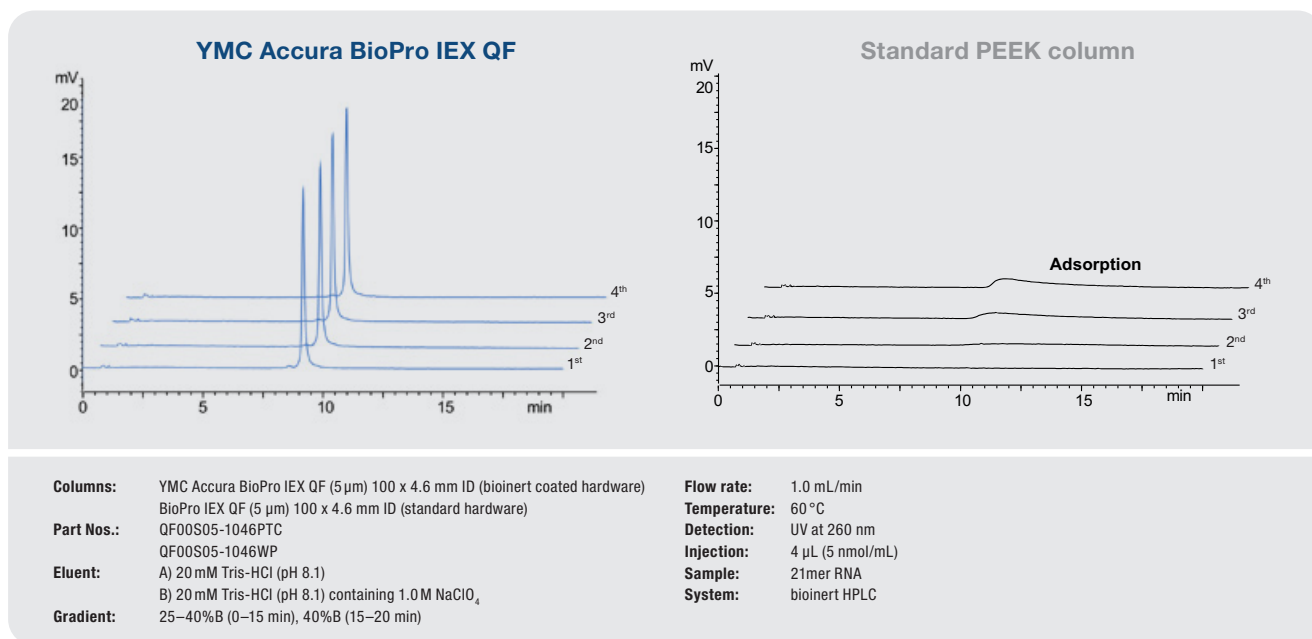

Oligonucleotides generally exhibit poor peak shape and therefore low recovery in AEX analysis, mainly due to adsorption onto the column hardware. YMC Accura BioPro IEX columns provide high recovery and very good peak shapes from the first injection. This makes YMC Accura BioPro IEX QF columns ideal for the analysis of oligonucleotides with reproducible results. The columns show stable peak areas from the first injection, so that no preconditioning is required.

YMC ACCURA

Different column lengths for each separation purpose

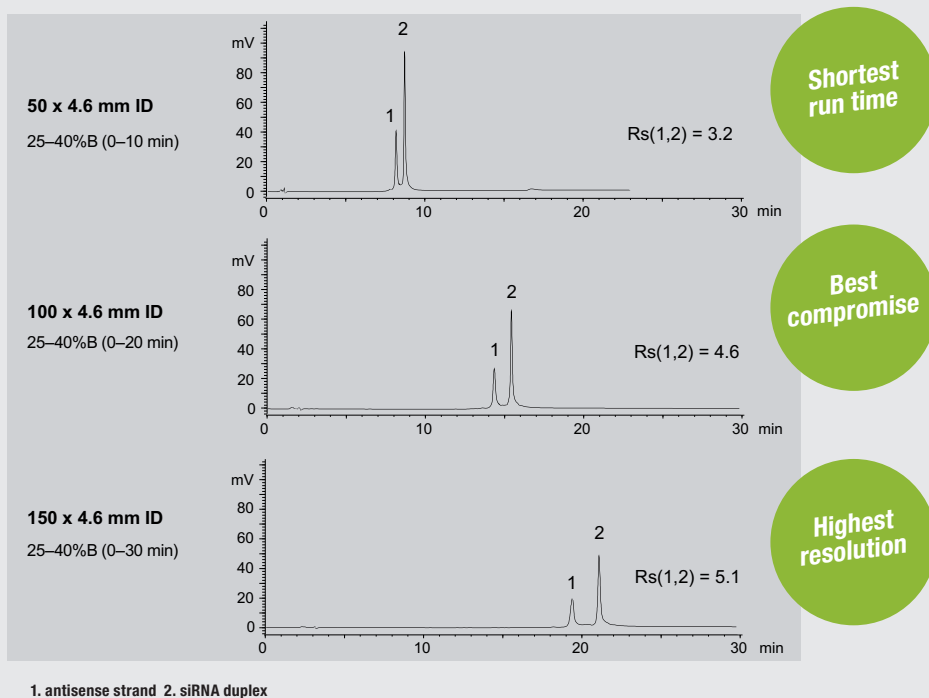

Columns: YMC Accura BioPro IEX QF (5 µm)

Part Nos.: QF00S05-0546PTC
QF00S05-1046PTC
QF00S05-1546PTC

Eluent: A) 20 mM Tris-HCl (pH 8.1)
B) 20 mM Tris-HCl (pH 8.1) containing 1.0 M NaClO₄

Flow rate: 0.5 mL/min

Temperature: 60 °C


Detection: UV at 260 nm

Injection: 8 µL (each 5 nmol/mL)

System: bioinert HPLC

AEX – STABILITY AND REPRODUCIBILITY


Excellent stability using temperatures up to 80 °C

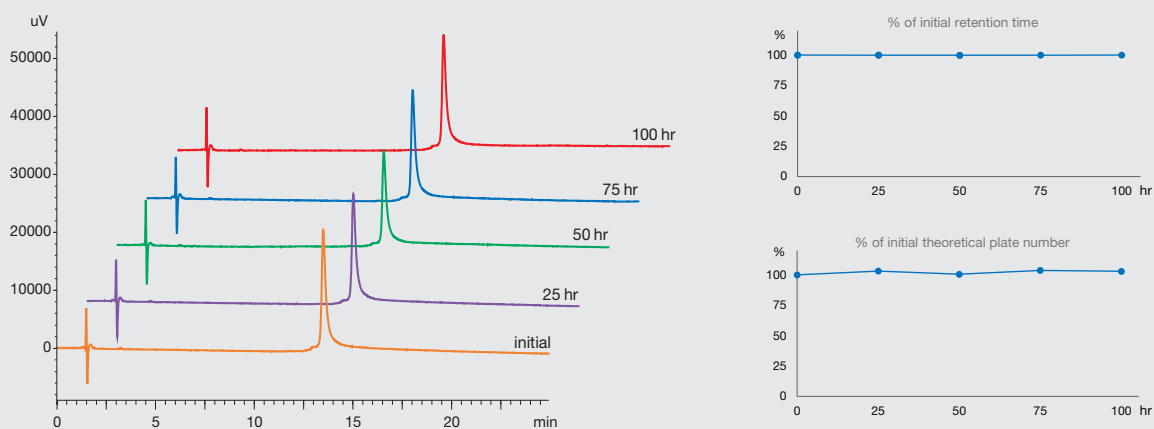

Column: YMC Accura BioPro IEX QF (5 µm) 100 x 4.6 mm ID
Part No.: QF00S05-1046PTC
Eluent: A) 20 mM Tris-HCl (pH 8.1)
 B) 20 mM Tris-HCl (pH 8.1) containing 1.0 M NaCl
Gradient: 40–90%B (0–25 min), 90%B (25–35 min), 40%B (35–60 min)

Flow rate: 0.5 mL/min
Temperature: 80 °C
Detection: UV at 260 nm
Injection: 4 µL (5 nmol/mL)
Sample: 21mer RNA




Excellent batch-to-batch reproducibility

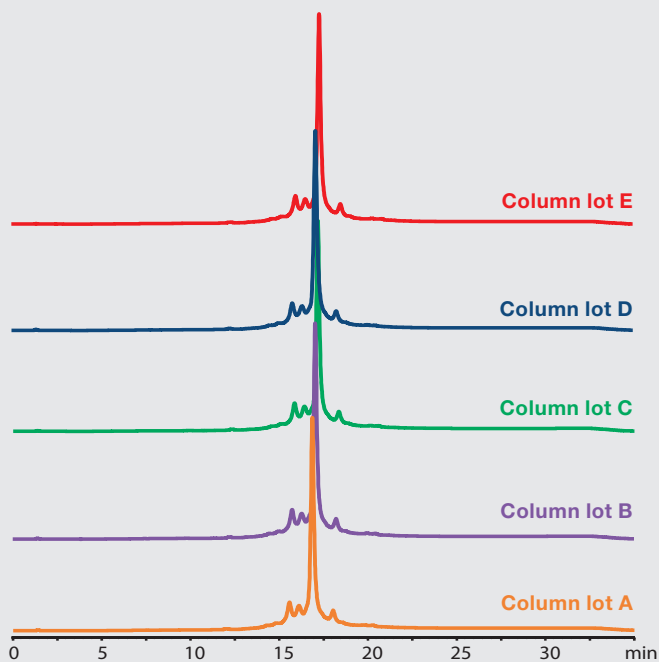

Column: BioPro IEX SF (5 µm) 100 x 4.6 mm ID
Part No.: SF00S05-1046WP
Eluent: A) 20 mM NaH₂PO₄ - Na₂HPO₄ (pH 6.5)
 B) 20 mM NaH₂PO₄ - Na₂HPO₄ (pH 6.5) containing 0.2 M NaCl
Gradient: 0–50%B (0.5–30 min)
Flow rate: 0.5 mL/min (180 cm/hr)
Temperature: 25 °C
Detection: UV at 215 nm
Injection: 10 µL
Sample: monoclonal antibody (IgG1)

YMC SEC

YMC COLUMN

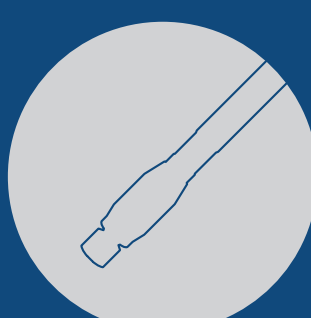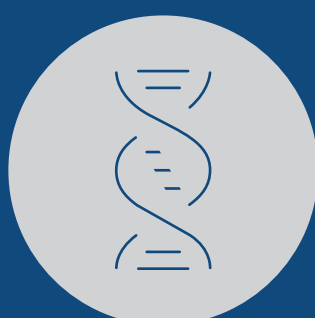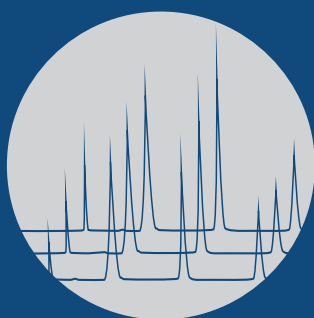


SEC – UHPLC / HPLC SELECTIVITIES

FEATURES

- Excellent reproducibility with minimal secondary interactions
- 2 µm for UHPLC
- Cost-effective

	YMC-Pack Diol-120	YMC-Pack Diol-200	YMC-Pack Diol-300
	For short oligonucleotides	For intermediate oligonucleotides	For longer oligonucleotides
Base particle	silica		
Particle size / µm	3, 5	2, 3, 5	2, 3, 5
Pore size / nm	12	20	30
Modification	dihydroxypropyl		
Temperature range	40 °C		
Pressure limit	2 µm: 45 MPa (6,525 psi); 3/5 µm: 20 MPa (3,000 psi)		


Column Selection Tool according to nucleotide length

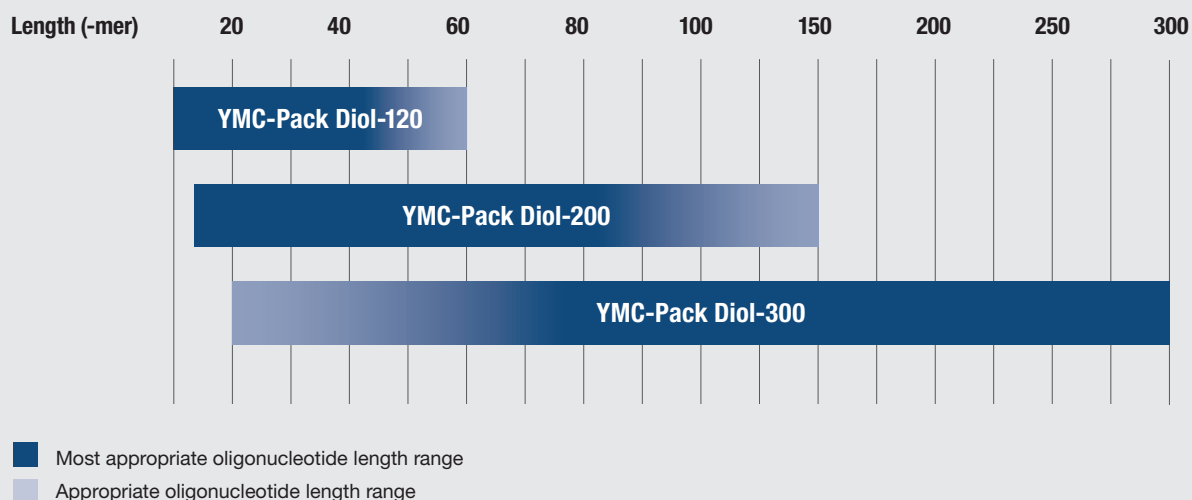


SEC – PORE SIZE SELECTION

SEC ANALYSIS USING YMC-PACK DIOL COLUMNS WITH DIFFERENT PORE SIZES

Analysis of ssDNA of 10–120mer length

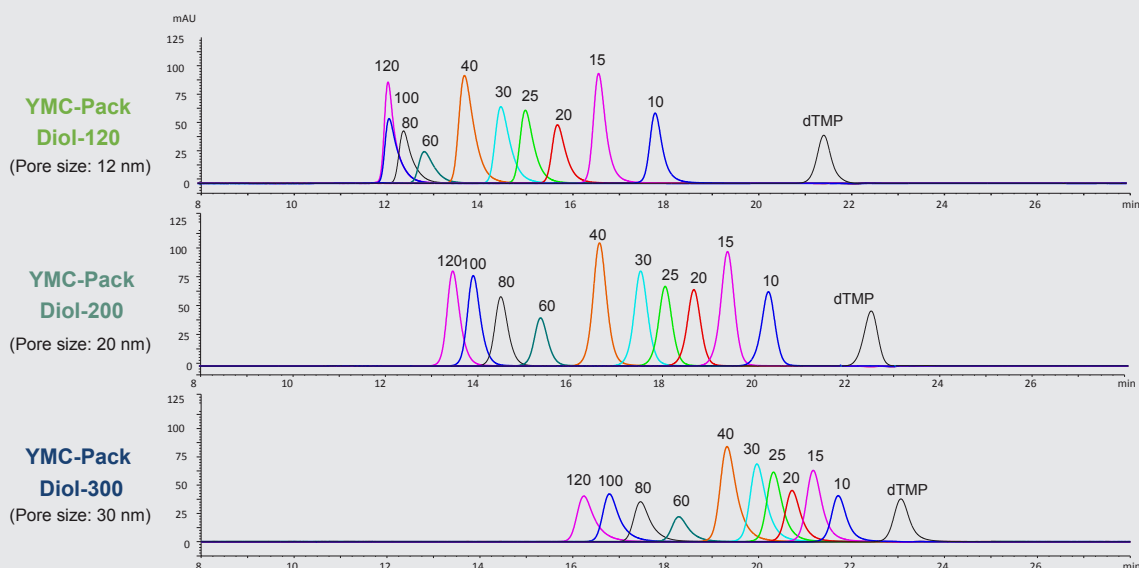

Analysis of dsDNA of 10–300mer length

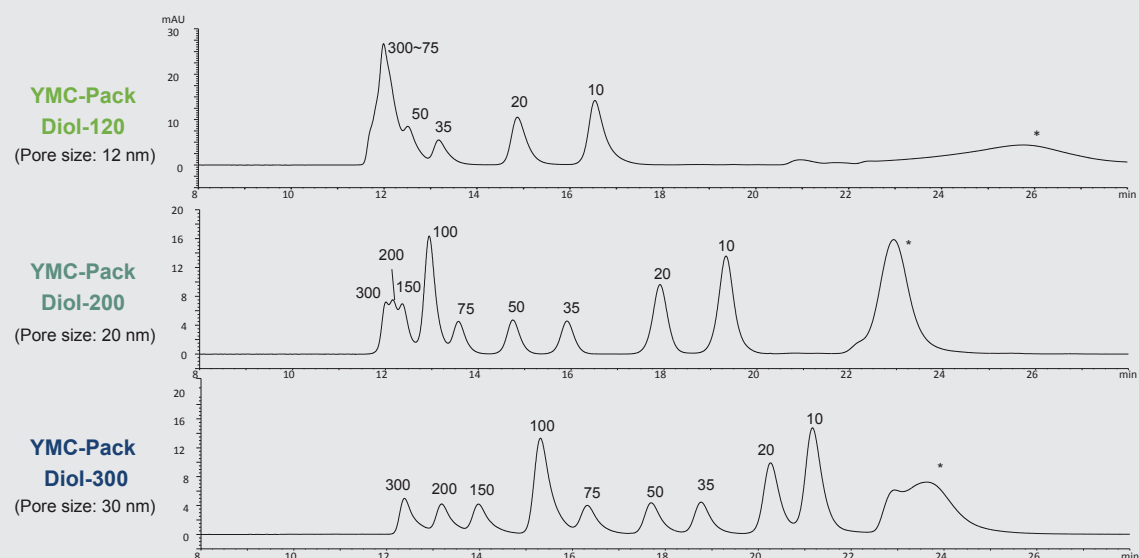

Columns: YMC-Pack Diol-120 (5 μ m, 12 nm)
300 x 4.6 mm ID
YMC-Pack Diol-200 (5 μ m, 20 nm)
300 x 4.6 mm ID
YMC-Pack Diol-300 (5 μ m, 30 nm)
300 x 4.6 mm ID

Part Nos.: DL12S05-3046WT
DL20S05-3046WT
DL30S05-3046WT
Eluent: 0.1 M KH_2PO_4 - K_2HPO_4 (pH 7.0) containing
0.2 M NaCl
Flow rate: 0.17 mL/min


Temperature: 25 $^{\circ}\text{C}$
Detection: UV at 260 nm
Injection: 1.0 μ L (each 5 nmol/mL)
Sample: ssDNA (10–120 mer + dTMP)
dsDNA (10–300 mer)

YMC-Pack Diol-120 is best suited for shorter oligonucleotides (10–40mer), whereas YMC-Pack Diol-200 shows the best resolution for oligonucleotides of medium size (30–80mer). Longer oligonucleotides of 60–120mer in length are separated most effectively by YMC-Pack Diol-300. Similar results are obtained when analysing dsDNA. Small oligonucleotides are separated with higher resolution when smaller pore sizes of 12 and 20 nm are used. Above a length of 50mer, oligonucleotides are unable to penetrate the small pores and elute at the same time. YMC-Pack Diol-200 can resolve oligonucleotides up to a size of 100mer. dsDNA of 150–300mer are only separated by YMC-Pack Diol-300 with the largest pore size of 30 nm. This column also shows the best resolution over a wide range of oligonucleotide lengths.

SEC – ELUTION VOLUMES

COMPARISON OF THE ELUTION VOLUME OF ssDNA AND dsDNA

Selected corresponding ssDNA and dsDNA pairs of the same number of bases are marked

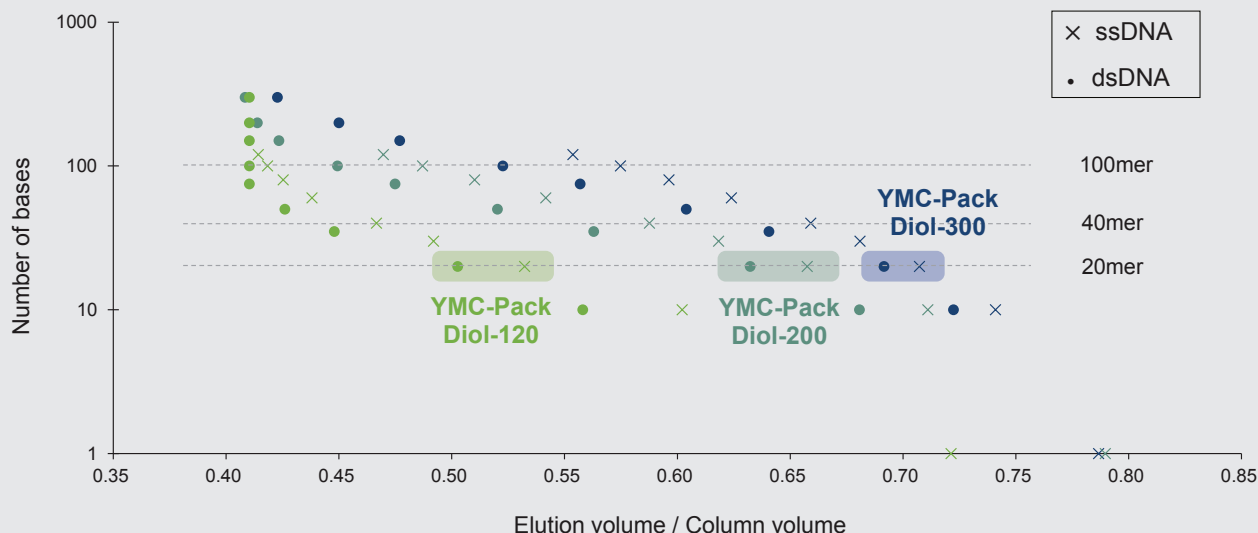

Columns: YMC-Pack Diol-120 (5 µm, 12 nm)
300 x 4.6 mm ID
YMC-Pack Diol-200 (5 µm, 20 nm)
300 x 4.6 mm ID
YMC-Pack Diol-300 (5 µm, 30 nm)
300 x 4.6 mm ID



Part Nos.: DL12S05-3046WT
DL20S05-3046WT
DL30S05-3046WT
Eluent: 0.1 M KH_2PO_4 - K_2HPO_4 (pH 7.0)
containing 0.2 M NaCl
Flow rate: 0.17 mL/min

Detection: UV at 260 nm
Temperature: 25 °C
Injection: 1.0 µL (each 5 nmol/mL)
Samples: ssDNA (10–120 mer + dTMP)
dsDNA (10–300 mer)

Although dsDNA has the same length as its single-stranded counterpart, the dsDNA elutes at lower elution volumes when separated by SEC. This behaviour is most probably due to the larger hydrodynamic radius of dsDNA compared to ssDNA, which results in faster diffusion through the stationary phase.

YMC HILIC

YMC COLUMN

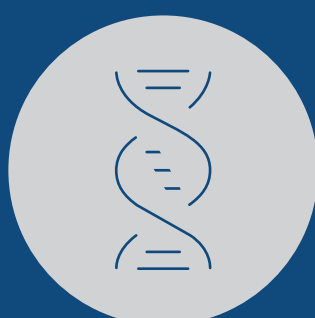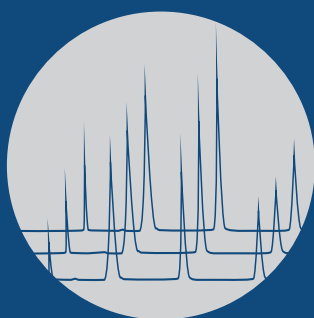


HILIC – UHPLC/HPLC SELECTIVITY

FEATURES


- pH- and temperature stable
- Superior reproducibility
- Bioinert coated YMC Accura column hardware

	Base particle	Modification	Particle Size/ μm	Pore Size/ nm	pH range	Temperature range	Column hardware
YMC-Triart Diol-HILIC	organic/inorganic hybrid silica	Diol (USP L20)	1.9, 3, 5	12	2–10	50 °C	bioinert coated YMC Accura


Pre-conditioning of a stainless-steel and a bioinert coated column with short DNA mixture

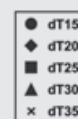

YMC-Triart Diol-HILIC (standard column)

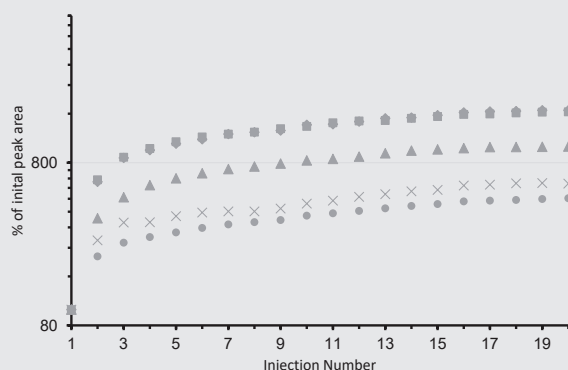

YMC Accura Triart Diol-HILIC

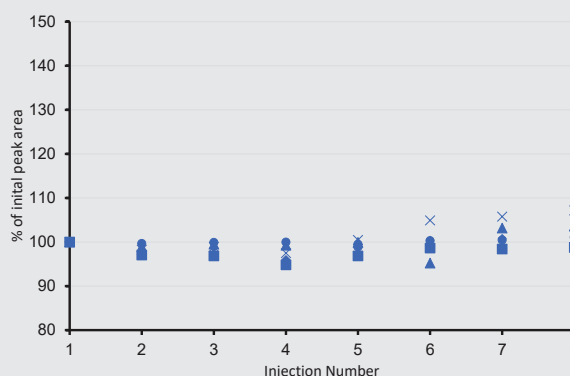

Columns: YMC-Triart Diol-HILIC (1.9 μm, 12 nm)
150 x 2.1 mm ID (standard hardware)
YMC Accura Triart Diol-HILIC (1.9 μm,
12 nm) 150 x 2.1 mm ID (bioinert hardware)

Part Nos.: TDH12SP9-15Q1WT
TDH12SP9-15Q1PTC

Eluent: A) 50 mM ammonium acetate (pH 6.9)
B) acetonitrile

Gradient: 75–45%B (0–30 min)

Flow rate: 0.3 mL/min

Temperature: 40 °C

Detection: UV at 260 nm


Injection: 2 μL

Sample: deoxythymidine oligonucleotides:
dT15-35 (2 μM)

System: bioinert system

Pre-conditioning is a typical procedure when working with stainless-steel columns. Using a bioinert column such as YMC Accura Triart usually achieves great performance from the first injection when working with an IP-RP phase. HILIC phases still need some pre-conditioning when a bioinert column is used; however, the number of injections is remarkably reduced. While 20 injections are necessary for the stainless-steel column, the YMC Accura column is already conditioned after 8 injections, with very little difference (less than 10 %) between initial and final peak areas.

Improved chromatographic results using bioinert coated
YMC Accura Triart column

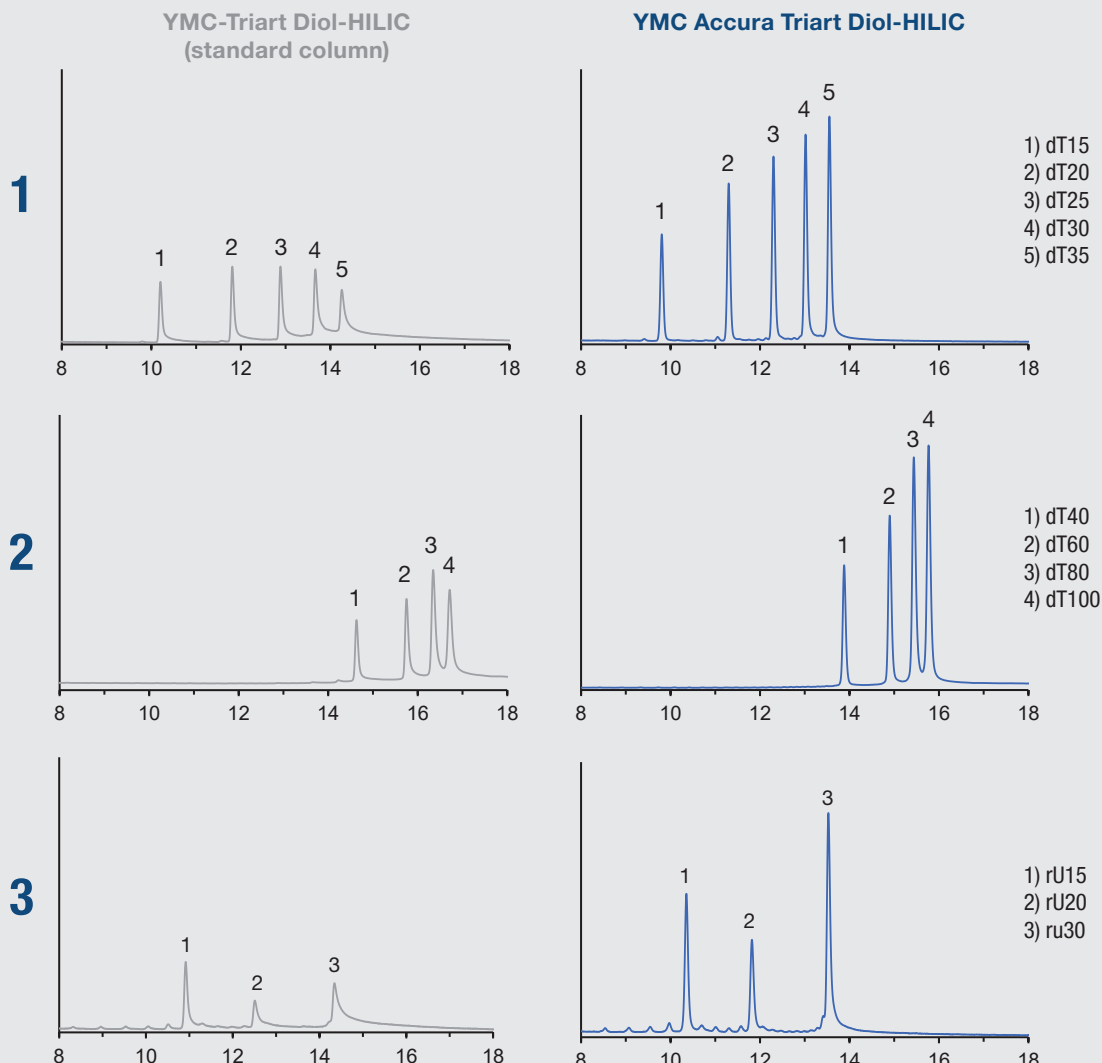

Columns: YMC-Triart Diol-HILIC (1.9 μ m, 12 nm)
150 x 2.1 mm ID (standard hardware)
YMC Accura Triart Diol-HILIC (1.9 μ m,
12 nm) 150 x 2.1 mm ID (bioinert hardware)
Part Nos.: TDH12SP9-15Q1WT
TDH12SP9-15Q1PTC

Eluent: A) 50 mM ammonium acetate (pH 6.9)
B) acetonitrile
Gradient: 75–45%B (0–30 min)
Flow rate: 0.3 mL/min
Temperature: 40 °C
Detection: UV at 260 nm


Injection: 2 μ L
Sample: deoxythymidine oligonucleotides:
dT15-35 (2 μ M) and dT40-100 (2 μ M)
RNA oligonucleotides: rU15-30 (2 μ M)
System: bioinert system

By courtesy of Honorine Lardeux, Institute of Pharmaceutical Sciences of Western Switzerland (University of Geneva), Geneva, Switzerland.
Reference: H. Lardeux, A. Goyon, K. Zhang, J.M. Nguyen, M.A. Lauber, D. Guilleme, V. D'Atri, The impact of low adsorption surfaces
for the analysis of DNA and RNA oligonucleotides, J. Chromatogr. A 1677 (2022) 463324.

After conditioning and analysing the short DNA oligonucleotide mixture of dT15-35, longer DNA oligonucleotides dT40-100 and short RNA oligonucleotides rU15-30 are analysed. Higher sensitivities, peak areas and less tailing are achieved using the bioinert YMC Accura Triart Diol-HILIC column. Non-specific adsorption does not vary according to length, even though the adsorption is usually higher for longer oligonucleotides in IP-RP.

HILIC EXPERT TIPS

Influence of sample solvent

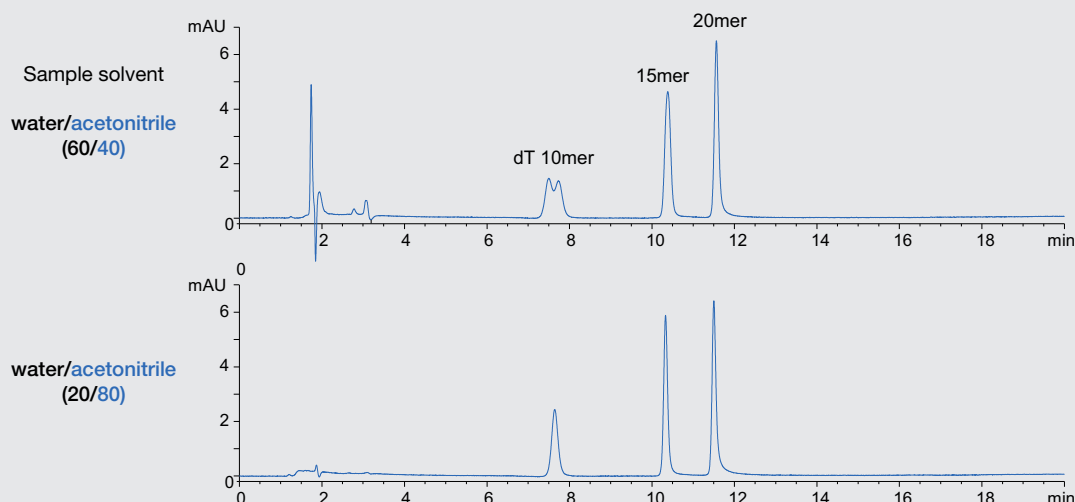


Column: YMC Accura Triart Diol-HILIC (3 μ m, 12 nm), 150 X 2.1 mm ID
Part No.: TDH12S03-15Q1PTC
Eluent: A) 50mM HCOONH₄ (pH 6.5)/acetonitrile (30/70)
 B) 50mM HCOONH₄ (pH 6.5)
Gradient: 0–43%B (0–20 min), 43%B (20–25 min)
 ((70–40% acetonitrile (0–20 min), 40% acetonitrile (20–25 min))

Flow rate: 0.21 mL/min
Temperature: 40 °CC
Sample: Oligo DNA (poly(dT) 10, 15, 20mer)
Injection: 2 μ L (each 2.0 nmol/mL)

The influence of the sample solvent on the peak shape is significant. The organic composition of the sample solvent must be equal to or higher than the initial gradient composition. A higher water content in the sample solvent leads to massive peak deformation.

Influence of mobile phase pH

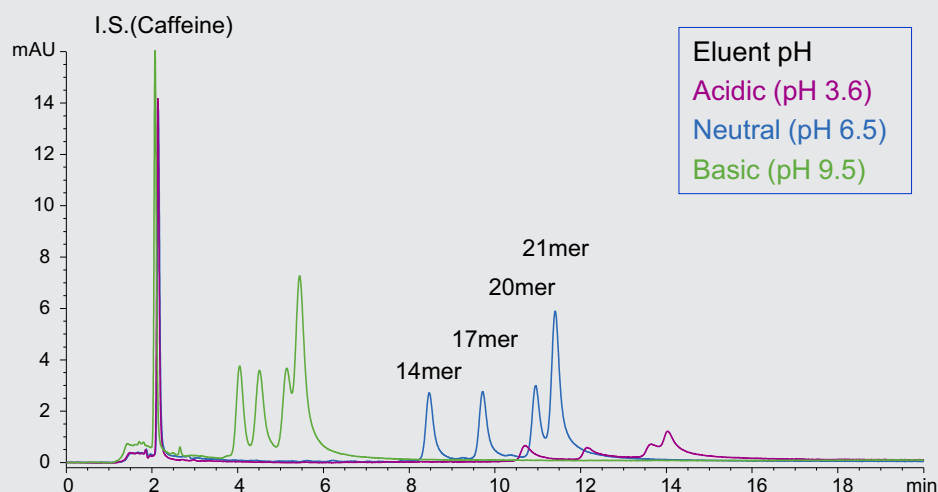



Column: YMC Accura Triart Diol-HILIC (3 μ m, 12 nm), 150 x 2.1 mm ID
Part No.: TDH12S03-15Q1PTC
Eluent: A) 50 mM ammonium formate buffer/acetonitrile (30/70)
 B) 50 mM ammonium formate buffer
Gradient: 7.1–21.4%B (0–20 min), 21.4%B (20–25 min)
 (65–55% acetonitrile (0–20 min), 55% acetonitrile (20–25 min))

Flow rate: 0.21 mL/min
Temperature: 40 °C
Injection: 1 μ L (1.8 nmol/mL-3.6 nmol/mL)
Sample: All PO RNA (14, 17, 20, 21mer) + I.S.(Caffeine)

The pH of the mobile phase has a massive effect on the recovery and retention of oligonucleotides. Acidic pH results in higher retention, but at the cost of drastically reduced recovery. A neutral-to-basic pH is recommended for the mobile phase. When using a basic pH, a shorter retention time is observed, but also the highest recovery.

YMC ORDERING INFORMATION

YMC COLUMN

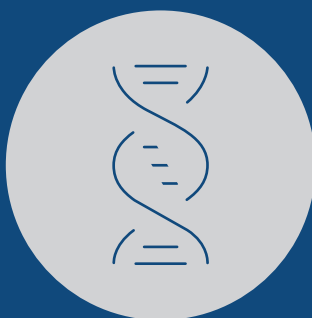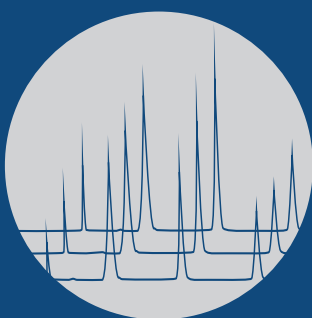

IP-RP – ORDERING INFORMATION

1.9 µm bioinert coated UHPLC columns (max. pressure 100 MPa)

Phase	Column ID [mm]	Column length [mm]					Guard cartridges* with 5 mm length
		20	30	50	100	150	pack of 3 (EMEA) 1 piece (outside EMEA)
YMC Accura Triart C18	2.1	TA12SP9-02Q1PTC	TA12SP9-03Q1PTC	TA12SP9-05Q1PTC	TA12SP9-10Q1PTC	TA12SP9-15Q1PTC	TA12SP9-E5Q1GCC
YMC Accura Triart Bio C18	2.1	TA30SP9-02Q1PTC	TA30SP9-03Q1PTC	TA30SP9-05Q1PTC	TA30SP9-10Q1PTC	TA30SP9-15Q1PTC	TA30SP9-E5Q1GCC
YMC Accura Triart C8	2.1	TO12SP9-02Q1PTC	TO12SP9-03Q1PTC	TO12SP9-05Q1PTC	TO12SP9-10Q1PTC	TO12SP9-15Q1PTC	TO12SP9-E5Q1GCC
YMC Accura Triart Bio C4	2.1	TB30SP9-02Q1PTC	TB30SP9-03Q1PTC	TB30SP9-05Q1PTC	TB30SP9-10Q1PTC	TB30SP9-15Q1PTC	TB30SP9-E5Q1GCC

3 µm bioinert coated HPLC columns (max. pressure 45 MPa)

Phase	Column ID [mm]	Column length [mm]			Guard cartridges* with 5 mm length
		50	100	150	pack of 3 (EMEA) 1 piece (outside EMEA)
YMC Accura Triart C18	2.1	TA12S03-05Q1PTC	TA12S03-10Q1PTC	TA12S03-15Q1PTC	TA12S03-E5Q1GCC
	4.6	TA12S03-0546PTC	TA12S03-1046PTC	TA12S03-1546PTC	TA12S03-E546GCC
YMC Accura Triart Bio C18	2.1	TA30S03-05Q1PTC	TA30S03-10Q1PTC	TA30S03-15Q1PTC	TA30S03-E5Q1GCC
	4.6	TA30S03-0546PTC	TA30S03-1046PTC	TA30S03-1546PTC	TA30S03-E546GCC
YMC Accura Triart C8	2.1	TO12S03-05Q1PTC	TO12S03-10Q1PTC	TO12S03-15Q1PTC	TO12S03-E5Q1GCC
	4.6	TO12S03-0546PTC	TO12S03-1046PTC	TO12S03-1546PTC	TO12S03-E546GCC
YMC Accura Triart Bio C4	2.1	TB30S03-05Q1PTC	TB30S03-10Q1PTC	TB30S03-15Q1PTC	TB30S03-E5Q1GCC
	4.6	TB30S03-0546PTC	TB30S03-1046PTC	TB30S03-1546PTC	TB30S03-E546GCC

*Guard cartridge holder required, part no. XPCHUHP

3 µm HPLC columns (max. pressure 20-30 MPa)

Phase	Column ID [mm]	Column length [mm]					Guard cartridges** with 10 mm length
		50	75	100	150	250	pack of 5
Hydrosphere C18	2.1	HS12S03-05Q1WT	HS12S03-L5Q1WT	HS12S03-10Q1WT	HS12S03-15Q1WT	HS12S03-25Q1WT	HS12S03-01Q1GC
	3.0	HS12S03-0503WT	HS12S03-L503WT	HS12S03-1003WT	HS12S03-1503WT	HS12S03-2503WT	HS12S03-0103GC
	4.6	HS12S03-0546WT	HS12S03-L546WT	HS12S03-1046WT	HS12S03-1546WT	HS12S03-2546WT	HS12S03-0104GC

5 µm HPLC columns (max. pressure 20–25 MPa)

Phase	Column ID [mm]	Column length [mm]					Guard cartridges** with 10 mm length
		50	75	100	150	250	pack of 5
Hydrosphere C18	2.1	HS12S05-05Q1WT	HS12S05-L5Q1WT	HS12S05-10Q1WT	HS12S05-15Q1WT	HS12S05-25Q1WT	HS12S05-01Q1GC
	3.0	HS12S05-0503WT	HS12S05-L503WT	HS12S05-1003WT	HS12S05-1503WT	HS12S05-2503WT	HS12S05-0103GC
	4.6	HS12S05-0546WT	HS12S05-L546WT	HS12S05-1046WT	HS12S05-1546WT	HS12S05-2546WT	HS12S05-0104GC

**Guard cartridge holder required, part no. XPGCH-Q1 (for EMEA)/XPGCHP1 (outside EMEA)

IP-RP – ORDERING INFORMATION

5 µm bioinert coated HPLC columns (max. pressure 10–45 MPa)

Phase	Column ID [mm]	Column length [mm]			Guard cartridges* with 5 mm length
		50	100	150	pack of 3 (EMEA) 1 piece (outside EMEA)
YMC Accura Triart C18	2.1	TA12S05-05Q1PTC	TA12S05-10Q1PTC	TA12S05-15Q1PTC	TA12S05-E5Q1GCC
	4.6	TA12S05-0546PTC	TA12S05-1046PTC	TA12S05-1546PTC	TA12S05-E546GCC
	10	–	TA12S05-1010PTC	TA12S05-1510PTC	–
YMC Accura Triart Bio C18	2.1	TA30S05-05Q1PTC	TA30S05-10Q1PTC	TA30S05-15Q1PTC	TA30S05-E5Q1GCC
	4.6	TA30S05-0546PTC	TA30S05-1046PTC	TA30S05-1546PTC	TA30S05-E546GCC
	10	–	TA30S05-1010PTC	TA30S05-1510PTC	–
YMC Accura Triart C8	2.1	T012S05-05Q1PTC	T012S05-10Q1PTC	T012S05-15Q1PTC	T012S05-E5Q1GCC
	4.6	T012S05-0546PTC	T012S05-1046PTC	T012S05-1546PTC	T012S05-E546GCC
	10	–	T012S05-1010PTC	T012S05-1510PTC	–
YMC Accura Triart Bio C4	2.1	TB30S05-05Q1PTC	TB30S05-10Q1PTC	TB30S05-15Q1PTC	TB30S05-E5Q1GCC
	4.6	TB30S05-0546PTC	TB30S05-1046PTC	TB30S05-1546PTC	TB30S05-E546GCC
	10	–	TB30S05-1010PTC	TB30S05-1510PTC	–

*Guard cartridge holder required, part no. XPCUHP

5 µm YMC-Actus high-throughput (semi)preparative columns (max. pressure 20–30 MPa)

Phase	Column ID [mm]	Column length [mm]					Guard cartridge** with 10 mm length
		50	75	100	150	250	pack of 2
YMC-Triart C18	20	TA12S05-0520WX	TA12S05-L520WX	TA12S05-1020WX	TA12S05-1520WX	TA12S05-2520WX	TA12S05-0120CCN
	30	TA12S05-0530WX	TA12S05-L530WX	TA12S05-1030WX	TA12S05-1530WX	TA12S05-2530WX	TA12S05-0130CCN
	50	TA12S05-0553DX	–	TA12S05-1053DX	TA12S05-1553DX	TA12S05-2553DX	TA12S05-0553DXG***
YMC-Triart Bio C18	20	TA30S05-0520WX	TA30S05-L520WX	TA30S05-1020WX	TA30S05-1520WX	TA30S05-2520WX	TA30S05-0120CCN
	30	TA30S05-0530WX	TA30S05-L530WX	TA30S05-1030WX	TA30S05-1530WX	TA30S05-2530WX	TA30S05-0130CCN
	50	TA30S05-0553DX	–	TA30S05-1053DX	TA30S05-1553DX	TA30S05-2553DX	TA30S05-0553DXG***
YMC-Triart C8	20	T012S05-0520WX	T012S05-L520WX	T012S05-1020WX	T012S05-1520WX	T012S05-2520WX	T012S05-0120CCN
	30	T012S05-0530WX	T012S05-L530WX	T012S05-1030WX	T012S05-1530WX	T012S05-2530WX	T012S05-0130CCN
	50	T012S05-0553DX	–	T012S05-1053DX	T012S05-1553DX	T012S05-2553DX	T012S05-0553DXG***
YMC-Triart Bio C4	20	TB30S05-0520WX	TB30S05-L520WX	TB30S05-1020WX	TB30S05-1520WX	TB30S05-2520WX	TB30S05-0120CCN
	30	TB30S05-0530WX	TB30S05-L530WX	TB30S05-1030WX	TB30S05-1530WX	TB30S05-2530WX	TB30S05-0130CCN
	50	TB30S05-0553DX	–	TB30S05-1053DX	TB30S05-1553DX	TB30S05-2553DX	TB30S05-0553DXG***
Hydrosphere C18	20	HS12S05-0520WX	HS12S05-L520WX	HS12S05-1020WX	HS12S05-1520WX	HS12S05-2520WX	HS12S05-0120CCN
	30	HS12S05-0530WX	HS12S05-L530WX	HS12S05-1030WX	HS12S05-1530WX	HS12S05-2530WX	HS12S05-0130CCN

**Guard cartridge holder required, part no. XPGHFSP20ID (20 mm ID) XPGHFSP30ID (30 mm ID)

***no holder required for 50 mm

Bioinert
semi-preparative
columns are
available on
request.

AEX – ORDERING INFORMATION

3 µm bioinert coated HPLC columns (max. pressure 15–30 MPa)

Phase	Column ID [mm]	Column length [mm]					Guard cartridges* with 5 mm length
		20	30	50	100	150	pack of 3 (EMEA) 1 piece (outside EMEA)
YMC Accura BioPro IEX QF	2.1 4.6	QF00S03-02Q1PTC –	QF00S03-03Q1PTC –	QF00S03-05Q1PTC QF00S03-0546PTC	QF00S03-10Q1PTC QF00S03-1046PTC	QF00S03-15Q1PTC QF00S03-1546PTC	QF00S03-E5Q1GCC QF00S03-E546GCC

5 µm bioinert coated HPLC columns (max. pressure 10–30 MPa)

Phase	Column ID [mm]	Column length [mm]				Guard cartridges* with 5 mm length
		50	100	150	250	pack of 3 (EMEA) 1 piece (outside)
YMC Accura BioPro IEX QF	2.1 4.6	QF00S05-05Q1PTC QF00S05-0546PTC	QF00S05-10Q1PTC QF00S05-1046PTC	QF00S05-15Q1PTC QF00S05-1546PTC	– QF00S05-2546PTC	QF00S05-E5Q1GCC QF00S05-E546GCC

*Guard cartridge holder required, part no. XPCHUHP

6 µm non-porous (bioinert coated) semiprep. columns (max. pressure 3-9 MPa)

Phase	Column ID [mm]	Column length [100 mm]	
		stainless steel	bioinert coated stainless steel
(YMC Accura) BioPro IEX QF	10 20 30	QF00S06-1010WT QF00S06-1020WT QF00S06-1030WT	QF00S06-1010PTC QF00S06-1020PTC QF00S06-1030PTC

SEC – ORDERING INFORMATION

2 µm UHPLC columns (max. pressure 45 MPa)

Phase	Column ID [mm]	Column length [mm]		Guard cartridges* with 10 mm length
		150	300	pack of 5
YMC-Pack Diol-200	4.6	DL20S02-1546WT	DL20S02-3046WT	DL20S02-0104GC
YMC-Pack Diol-300	4.6	DL30S02-1546WT	DL30S02-3046WT	DL30S02-0104GC

*Guard cartridge holder required, part no. XPGCH-Q1 (for EMEA)/XPGCHP1 (outside EMEA)

3 µm HPLC columns (max. pressure 20 MPa)

Phase	Column ID [mm]	Column length [mm]			Guard cartridges* with 10/30 mm length
		150	250	300	pack of 5
YMC-Pack Diol-120	4.6	DL12S03-1546WT	DL12S03-2546WT	DL12S03-3046WT	DL12S03-0104GC
	6.0	–	–	DL12S03-3006WT	–
	8.0	DL12S03-1508WT	–	DL12S03-3008WT	DL12S03-0308WTG**
YMC-Pack Diol-200	4.6	DL20S03-1546WT	DL20S03-2546WT	DL20S03-3046WT	DL20S03-0104GC
	6.0	–	–	DL20S03-3006WT	–
	8.0	DL20S03-1508WT	–	DL20S03-3008WT	DL20S03-0308WTG**
YMC-Pack Diol-300	4.6	DL30S03-1546WT	DL30S03-2546WT	DL30S03-3046WT	DL30S03-0104GC
	6.0	–	–	DL30S03-3006WT	–
	8.0	DL30S03-1508WT	–	DL30S03-3008WT	DL30S03-0308WTG**

*Guard cartridge holder required, part no. XPGCH-Q1 (for EMEA)/XPGCHP1 (outside EMEA)

**no holder required for 30 x 8 mm ID guard columns (1 piece)
recommended column coupler part no. XRCP1602

5 µm HPLC columns (max. pressure 20 MPa)

Phase	Column ID [mm]	Column length [mm]		Guard cartridges* with 10/30 mm length
		250	300	pack of 5
YMC-Pack Diol-120	4.6	DL12S05-2546WT	DL12S05-3046WT	DL12S05-0104GC
	6.0	DL12S05-2506WT	DL12S05-3006WT	–
	8.0	–	DL12S05-3008WT	DL12S05-0308WTG**
	10.0	DL12S05-2510WT	DL12S05-3010WT	DL12S05-0310WTG**
YMC-Pack Diol-200	4.6	DL20S05-2546WT	DL20S05-3046WT	DL20S05-0104GC
	6.0	DL20S05-2506WT	DL20S05-3006WT	–
	8.0	–	DL20S05-3008WT	DL20S05-0308WTG**
	10.0	DL20S05-2510WT	DL20S05-3010WT	DL20S05-0310WTG**
YMC-Pack Diol-300	4.6	DL30S05-2546WT	DL30S05-3046WT	DL30S05-0104GC
	6.0	DL30S05-2506WT	DL30S05-3006WT	–
	8.0	–	DL30S05-3008WT	DL30S05-0308WTG**
	10.0	DL30S05-2510WT	DL30S05-3010WT	DL30S05-0310WTG**

*Guard cartridge holder required, part no. XPGCH-Q1 (for EMEA)/XPGCHP1 (outside EMEA)

**no holder required for 30 x 8 mm ID guard columns (1 piece)
recommended column coupler part no. XRCP1602 (for 8 mm ID) and XRCP1605 (for 10 mm ID)

**Bioinert
SEC columns
are available
on request.**

HILIC – ORDERING INFORMATION

1.9 µm bioinert coated UHPLC columns (max. pressure 100 MPa)

Phase	Column ID [mm]	Column length [mm]			Guard cartridges* with 5 mm length
		50	100	150	pack of 3 (EMEA) 1 piece (outside EMEA)
YMC Accura Triart Diol-HILIC	2.1	TDH12SP9-05Q1PTC	TDH12SP9-10Q1PTC	TDH12SP9-15Q1PTC	TDH12SP9-E5Q1GCC

3 µm bioinert coated HPLC columns (max. pressure 45 MPa)

Phase	Column ID [mm]	Column length [mm]			Guard cartridges* with 5 mm length
		50	100	150	pack of 3 (EMEA) 1 piece (outside EMEA)
YMC Accura Triart Diol-HILIC	2.1	TDH12S03-05Q1PTC	TDH12S03-10Q1PTC	TDH12S03-15Q1PTC	TDH12S03-E5Q1GCC
	4.6	TDH12S03-0546PTC	TDH12S03-1046PTC	TDH12S03-1546PTC	TDH12S03-E546GCC

5 µm bioinert coated HPLC columns (max. pressure 45 MPa)

Phase	Column ID [mm]	Column length [mm]			Guard cartridges* with 5 mm length
		50	100	150	pack of 3 (EMEA) 1 piece (outside EMEA)
YMC Accura Triart Diol-HILIC	2.1	TDH12S05-05Q1PTC	TDH12S05-10Q1PTC	TDH12S05-15Q1PTC	TDH12S05-E5Q1GCC
	4.6	TDH12S05-0546PTC	TDH12S05-1046PTC	TDH12S05-1546PTC	TDH12S05-E546GCC


*Guard cartridge holder required, part no. XPCHUHP

FURTHER DIMENSIONS AND GUARD CARTRIDGES
AVAILABLE IN REGULAR STAINLESS-STEEL HARDWARE.




The following brands, trademarks, or service marks are the property of the listed company and/or its subsidiaries.
Acquity, BEH, XBridge are trademarks of Waters Corp.
AdvanceBio is a trademark of Agilent Technologies Inc.
DNAPac is a trademark of Thermo Fisher Scientific Inc.
Every effort has been taken to ensure this list is accurate at the time of creating this brochure.

DOWNLOAD THE FOLLOWING DOCUMENTS

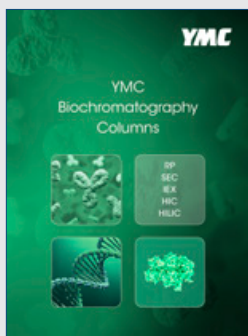


YMC Biochromatography Columns

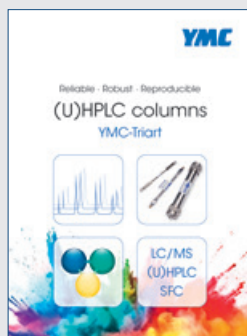

YMC-Triart



YMC BioPro IEX Resins

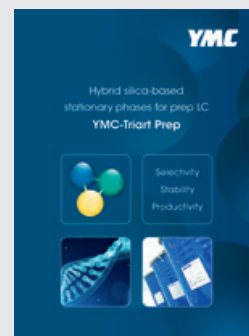


YMC-Triart Prep

YMC Brochures




Analysis and Purification of Oligonucleotides by AEX and IP-RP

YMC White Paper

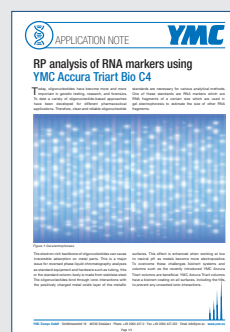


RP analysis of RNA markers using YMC Accura Triart Bio C4


YMC Application Notes

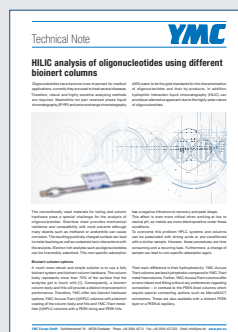

HILIC analysis of oligonucleotides using different bioinert columns

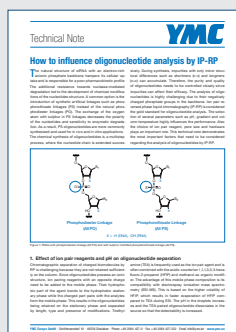

How to influence oligonucleotide analysis by IP-RP

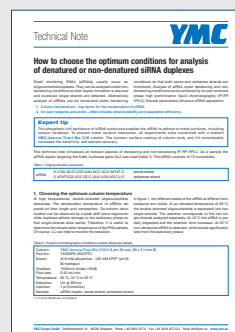

How to choose the optimum conditions for analysis of denatured or non-denatured siRNA duplexes

YMC Technical Notes



GET IN TOUCH
AND ORDER
NOW!

YMC COLUMN

Your local contact:



YMC CO., LTD.
www.ymc.co.jp

YMC Europe GmbH
www.ymc.eu

YMC America, Inc.
www.ymcamerica.com

YMC Schweiz GmbH
www.ymc-schweiz.ch

YMC India Pvt. Ltd.
www.ymcindia.com

YMC Biotechnology (Changzhou) Co., Ltd.

YMC Korea Co., Ltd.
www.ymckorea.com

YMC Taiwan Co., Ltd.
www.ymctaiwan.com